# Molecular Probes to Characterize HPLC Column Performance

Richard A. Henry<sup>1</sup>, Stephanie Schuster<sup>2</sup>, Conner McHale<sup>2</sup> and William Johnson<sup>2</sup>

<sup>1</sup>Independent Consultant, State College, PA 16801;

<sup>2</sup>Advanced Materials Technology Inc., Wilmington, DE



Presented at Pittcon 2019 Poster 1240-2

#### **Use of Chemical Probes in HPLC**


- "... the best overall characterization of a reversed-phase bonded phase is done by well-designed chromatographic tests." Uwe Neue, HPLC Columns, Wiley (1997)<sup>1.</sup> Probe mixtures have been developed by academic and industrial scientists since about 1990 in attempts to better understand HPLC column separations. Suppliers make use of probes to create columns that separate important classes of compounds. 1.Uwe D. Neue, HPLC Columns, Wiley-VCH (1997).
- User groups should consult probe studies published by researchers and column manufacturers extensively during HPLC method development. According to Neue, probes are the best way to identify (or eliminate) columns and mobile phases for selectivity toward target functional groups, and column manufacturers and academicians have best access to a variety of probe molecules. It is important to publish selectivity studies.
- Methanol is a particularly important mobile phase solvent. It should become more popular with users because column selectivity is often different and may be improved over acetonitrile in binary or ternary combinations with water.
- Superficially porous particle (SPP) columns facilitate high sample throughput. More columns and solvent systems can be evaluated to find optimum rather than adequate solutions. Lower pressure than sub-2 μm columns permits use of faster flow rates and higher viscosity methanol/water mobile phases. High efficiency and resolution minimizes the chance of peak overlap.

#### HALO® Reversed-Phase Columns for HPLC and UHPLC


	Packing Description	USP Code	Bonded Phase	Types of Interactions
	C18	L1	C18 (dimethyloctadecylsilane)	• Hydrophobic
l nc	C30	L62	C30 (Triacontyldimethylsilane)	• Hydrophobic
Increasing	С8	L7	C8 (dimethyloctylsilane)	• Hydrophobic
	AQ-C18	L1	proprietary	<ul> <li>Mainly hydrophobic</li> <li>Some dipole-dipole</li> </ul>
Polarity (s	Phenyl-Hexyl	L11	Phenyl-Hexyl (dimethylphenylhexylsilane)	• Hydrophobic • π - π
(solute	Biphenyl	L11	Biphenyl (dimethylbiphenylsilane)	• π – π • Hydrophobic
	RP-Amide	L60	C16 Amide	<ul><li>Hydrogen Bonding</li><li>Hydrophobic</li></ul>
dependent)	ES-CN	L10	ES-CN (di- <u>isopropylcyanopropylsilane</u> )	<ul><li>Dipole-dipole</li><li>Hydrophobic</li></ul>
ent)	PFP	L43	PFP (pentafluorophenylpropylsilane)	<ul> <li>Hydrophobic</li> <li>π - π</li> <li>Dipole-dipole</li> <li>Hydrogen bonding</li> </ul>


#### **Aromatic Compounds (with Pi electrons)**

\* <u>Pi electrons</u>- aromatic systems can be viewed as a large, planar surface with potential for polar interactions, such as Hydrogen Bond accepting ability. Delocalized electrons in the ring system interact with individual polar atoms and other aromatic networks to form strong donor-acceptor complexes that create unique chromatographic retention. Dissociation energies are similar in strength to H-Bonds.

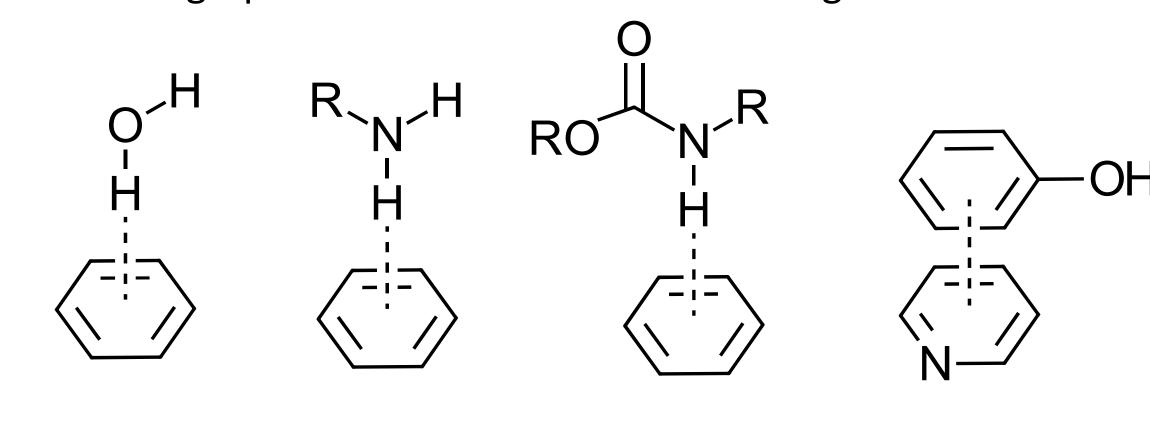

# Hydrophobic Subtraction Model<sup>2</sup>

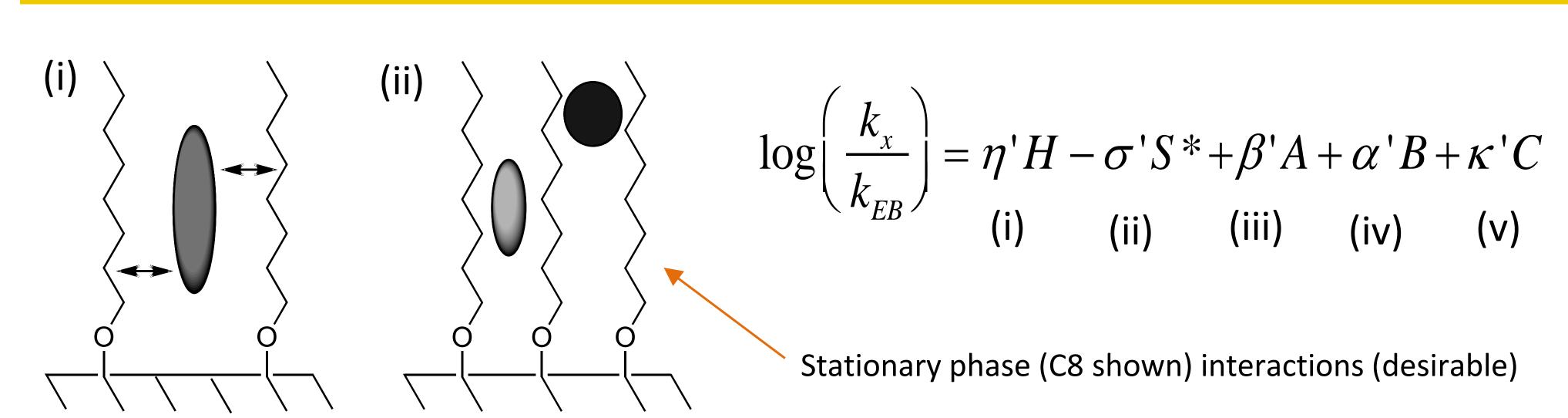

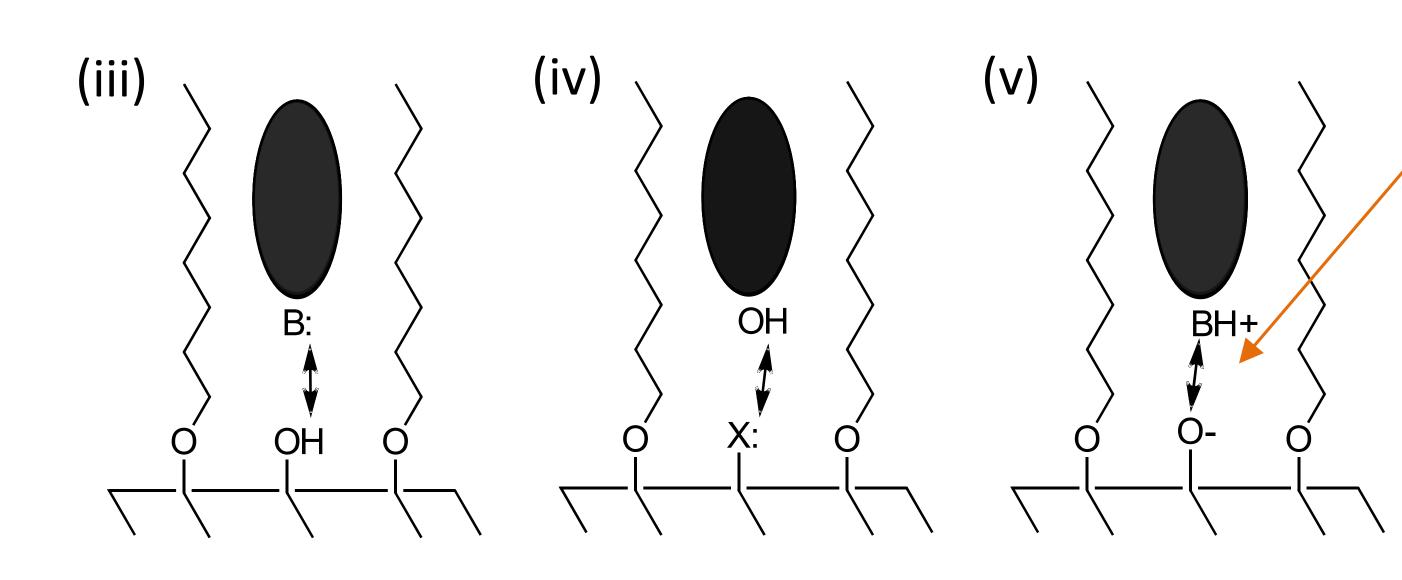

Silica substrate interactions

-Database available for 700 commercial columns, but equivalency comparisons are limited to MeCN mobile phases

2. Lloyd R. Snyder, Joseph J. Kirkland and John W. Dolan, *Introduction to Modern Liquid Chromatography*, John Wiley and Sons, Inc. (2010).

# Orthogonal HALO® Phases: Fs Values for HALO® Phases vs. HALO® C18


$$F_{S} = \sqrt{\left(w_{\mathbf{H}}(\mathbf{H}_{1} - \mathbf{H}_{2})\right)^{2} + \left(w_{\mathbf{S}^{*}}(\mathbf{S}^{*}_{1} - \mathbf{S}^{*}_{2})\right)^{2} + \left(w_{\mathbf{A}}(\mathbf{A}_{1} - \mathbf{A}_{2})\right)^{2} + \left(w_{\mathbf{B}}(\mathbf{B}_{1} - \mathbf{B}_{2})\right)^{2} + \left(w_{\mathbf{C}_{2.8}}(\mathbf{C}_{2.8_{1}} - \mathbf{C}_{2.8_{2}})\right)^{2}}$$

Fs	Name	Н	S*	A	В	C (pH 2.8)	C (pH 7.0)	ethylbenzene k
0	HALO C18	1.10	0.04	0	-0.05	0.05	0.04	6.1
12.07	HALO AQ-C18	1.00	-0.036	0.099	-0.048	0.156	0.864	6.7
10.04	HALO C8	0.91	0.02	-0.13	0	-0.01	0.18	4.3
17.43	HALO C30	0.938	-0.046	-0.14	0.023	0.17	0.35	4.5
52.83	HALO RP-Amide	0.85	0.08	-0.38	0.19	-0.41	0.31	4.6
17.35	HALO Phenyl- Hexyl	0.780	-0.09	-0.23	0	0.1	0.45	3.5
26.76	HALO Biphenyl	0.708	-0.183	-0.279	0.028	0.047	0.99	3.1
94.45	HALO PFP	0.702	-0.117	-0.073	-0.062	1.17	0.972	2.3
22.78	HALO ES-CN	0.566	-0.11	-0.344	0.021	0.126	1.15	1.88

- $F_s \le 3$ : columns are equivalent
- $3 < F_s < ^10$ : columns are likely equivalent for samples with a few
- $F_s \ge 35$ : columns are orthogonal

Data provided by Dwight Stoll group at Gustavus Adolphus College. www.hplccolumns.org

#### Impact of Phase Polarity for OH Probes

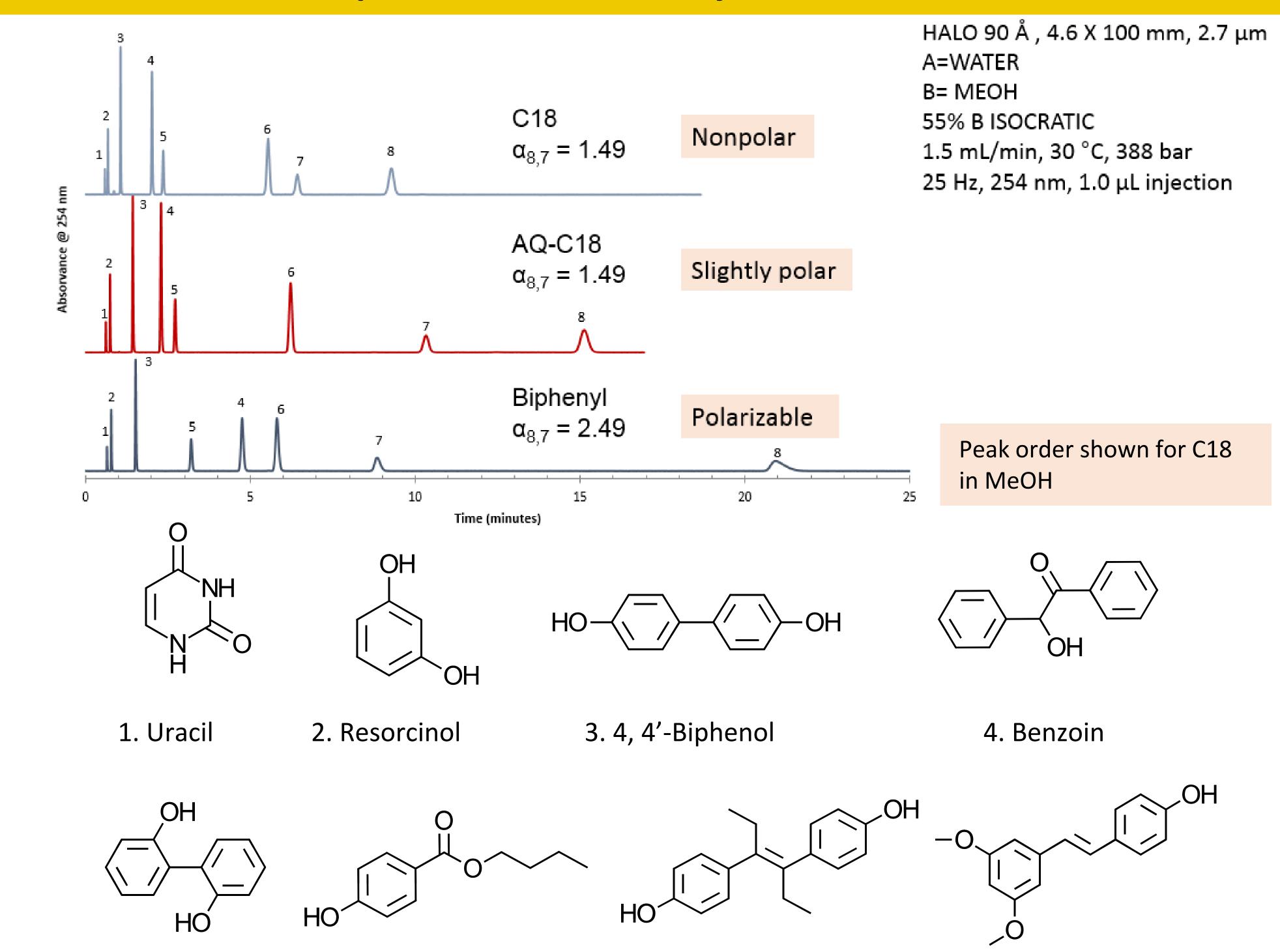

Retention and selectivity increases for solutes with strong OH dipoles although the stationary
phases above are not considered to have H-bonding ability; AQ-C18 has a small amount of dipole
character to increase wettability in pure water.

7. Diethylstilbesterol

Aromatic stationary phases, such as Biphenyl, show extra retention and even peak reversal for OH solutes because delocalized electrons are polarizable and more shape selective. Biphenyl is typically more retentive than Phenyl-Hexyl with different selectivity because it has two ring systems and no alkyl character.

#### **Enhanced Selectivity Values for OH Probe Mix in MeOH Mobile Phases**

HALO® C18	1.0	0
HALO® AQ-C18	0.970	17.3
HALO® C30	0.992	8.9
HALO® C8	0.880	34.6
HALO® Phenyl-Hexyl	0.971	17.0
HALO <sup>®</sup> Biphenyl	0.872	35.8
HALO® RP-Amide	0.836	40.4
DI	m.0	A 1 / 1
Phase in MeCN	R <sup>2</sup>	S Value
Phase in MeCN	R <sup>2</sup>	S Value
HALO® C18	1.0	S Value 0
HALO® C18	1.0	0
HALO® C18 HALO® AQ-C18	1.0 0.987	0 11.8
HALO® C18 HALO® AQ-C18 HALO® C30	1.0 0.987 0.995	0 11.8 7.1
HALO® C18 HALO® AQ-C18 HALO® C30 HALO® C8	1.0 0.987 0.995 0.979	0 11.8 7.1 14.5
HALO® C18 HALO® AQ-C18 HALO® C30 HALO® C8 HALO® Phenyl-Hexyl	1.0 0.987 0.995 0.979 0.988	0 11.8 7.1 14.5 11.0

Phase in MeOH

5. 2,2'-Biphenol 6. Butyl hydroxybenzoate


 $S = 100 V(1-R^2)$ 

Selectivity difference values by Neue are often enhanced in MeOH although the use of MeCN is more common because of lower operating pressure. One exception to this rule is often RP-Amide.


8. Pterostilbene

Using SPP particles, elevated temperatures or ternary MeCN/MeOH blends reduce operating pressure.

# Impact of Hydrophobicity Plus Strong Hydrogen Bonding

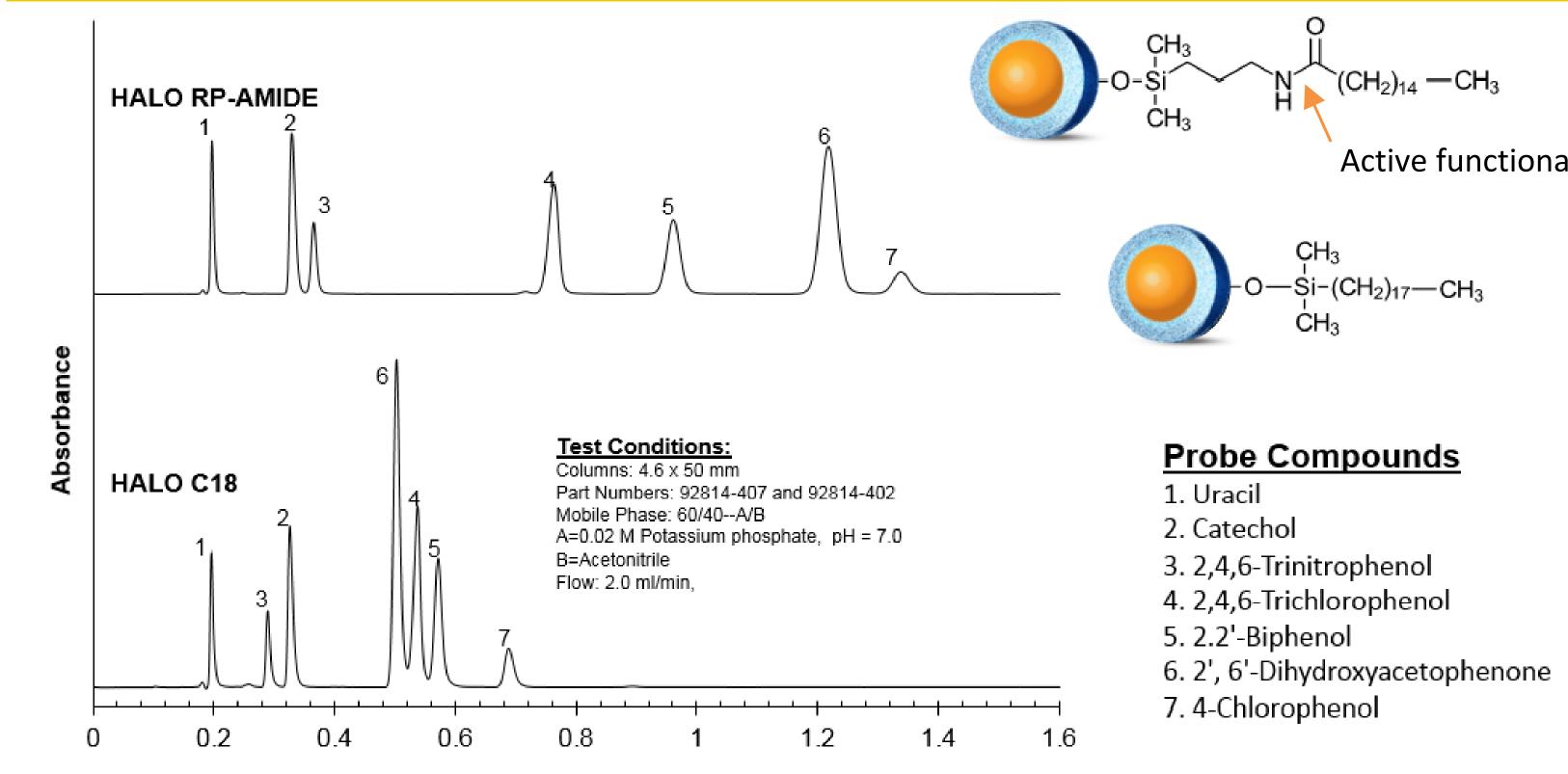

### **RP-Amide Shows Different Selectivity in MeOH for OH Probe**

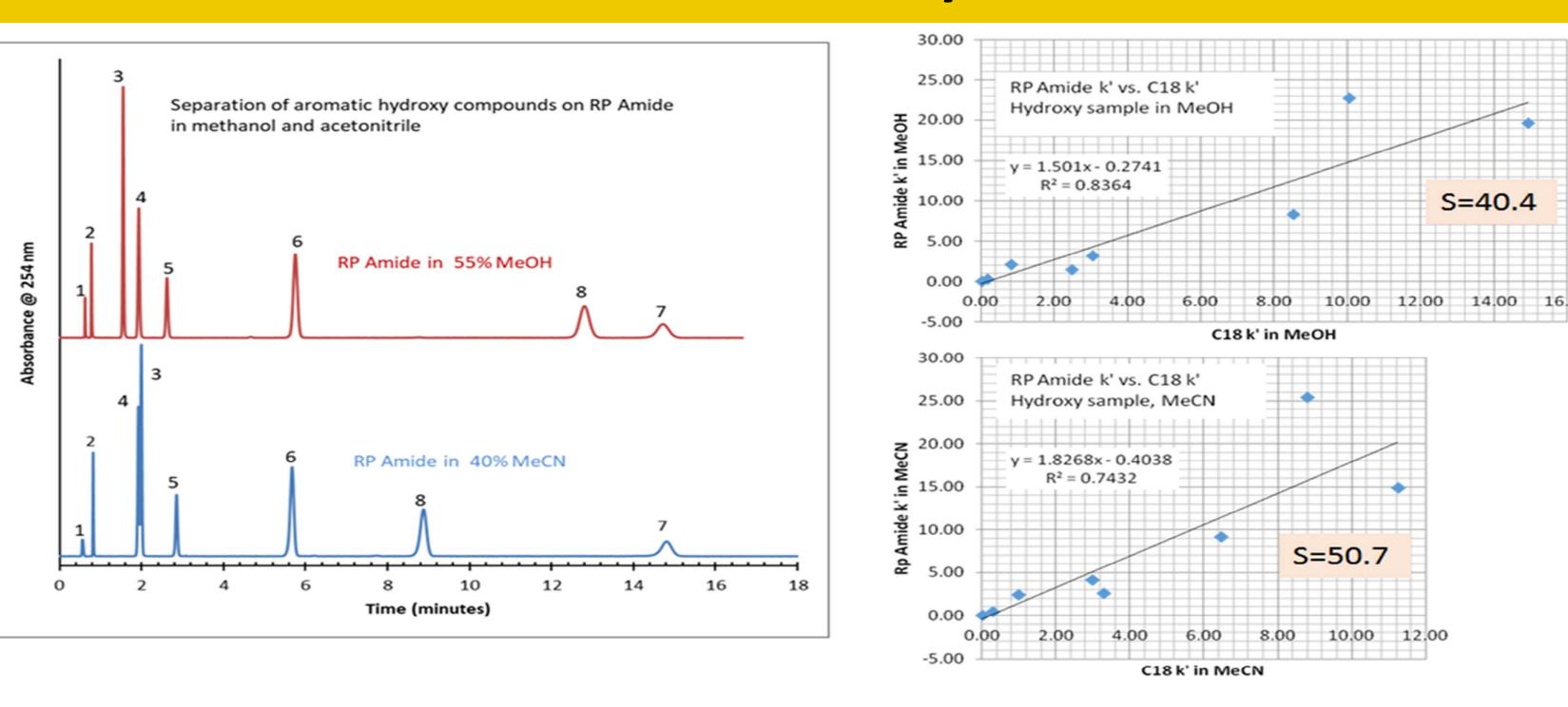

 While use of HALO® RP-Amide in MeCN mobile phase is more common, peak spacing and resolution may be improved in MeOH, making it preferred for this sample. Since H-bonding selectivity is different for RP-Amide in the two solvents, a ternary mobile phase may be ideal for difficult separations.

#### Probe Conclusions/ Acknowledgements

- Methanol is an important HPLC solvent for mobile phase use, but may be overlooked because
  acetonitrile works so well with C18 columns that are deeply entrenched in HPLC methods. In the
  interest of saving time, users are tempted to settle for adequate methods on C18 columns using
  acetonitrile mobile phases.
- Users should screen columns with sample components and the widest range of bonded phases that includes alternatives to hydrophobic interactions and not be limited to C18 and acetonitrile.
- Test probes are clearly the best way to develop a better understanding of how columns and mobile phases work, and skilled users should be more widely encouraged to publish results with probe compounds that are not proprietary.
- Rugged columns with SPP particles are ideal for broad early screening because their fundamental design guarantees fast separation results, moderate pressure, and high peak resolution. Time saved by screening with SPP columns allows study of more phases and variables such as methanol mobile phases. The lower operating pressure and high efficiency of SPP particles also makes SPP columns attractive for the final method.

-Acknowledgements to Mark Schure of Kroungold Consulting and to Advanced Materials Technology, Inc. for providing data and support during poster preparation.

-® HALO and Fused-Core are registered trademarks of Advanced Materials Technology, Inc

