

HALO® SMALL MOLECULE COLUMN CARE & USE SHEET



Description

HALO® columns are available in multiple superficially porous particle sizes, pore sizes, and stationary phases.

- Applicable to 1 to 10 mm ID columns

Operation Guidelines

- The direction of flow is marked on the column label.
- Water and all common organic solvents are compatible with HALO® columns.
- Running the column in the reversed flow direction is acceptable for all particle sizes and dimensions except for 2 µm columns.
- 2 µm HALO® columns are rated to 1000 bar.
- 2.7 and 5 µm HALO® columns are rated to 600 bar.

Small Molecule HALO® Phases					
Phase	Shipping Solvent	Storage Solvent	low pH/Max. Temp.	high pH/Max.Temp.	100% Aqueous Compatible
AQ-C18	ACN	ACN	2/60 °C	9/40 °C	√
C8	ACN/water	ACN	2/60 °C	9/40 °C	
C18	ACN/water	ACN	2/60 °C	9/40 °C	
LPH-C18	ACN/Water	ACN	1/90 °C	8/40 °C	
PCS	ACN	ACN	2/40 °C	7/60 °C	
C30	ACN/water	ACN	2/60 °C	9/40 °C	√
Phenyl-Hexyl	ACN/water	ACN	2/60 °C	9/40 °C	√
Biphenyl	Methanol	Methanol	2/60 °C	9/40 °C	√
PFP	ACN	ACN	2/60 °C	8/40 °C	√
ES-CN	ACN	ACN	1/80 °C	8/40 °C	√
RP-Amide	ACN/water	ACN	2/60 °C	9/40 °C	√
HILIC	ACN	ACN	1/60 °C	8/40 °C	√
Penta-HILIC	90/10 ACN/water	ACN	2/60 °C	9/40 °C	√

*All small molecule phases have 90 Å pore size except C30 which has a pore size of 160 Å

Column Care & Storage

To maximize column life, ensure that samples and mobile phases are particle-free. The use of an in-line filter with 0.5-micron porosity between the pump and sample injector is highly recommended. Before storing the column, the end-fittings should be tightly sealed with the end-plugs that came with the column to prevent the packing from drying. For storage longer than 3 days, rinse the column with at least 20 column volumes of 50/50 ACN/water (if using buffer) then with 100% of the storage solvent. For column cleaning procedures, please see the HALO® Column Cleaning Procedures and Best Practices document by using the QR code.

Safety

- HPLC columns are for laboratory use only. Not for drug, household, or other use.
- Users of HPLC columns should be aware of the toxicity or flammability of the mobile phases chosen for use with the columns. Precautions should be taken to avoid contact and leaks.
- HPLC columns should be used in well-ventilated environments to minimize concentration of solvent fumes.

Ordering Information & Technical Support

For ordering information or for technical support on this product, please contact us at support@advanced-materials-tech.com or halocolumns.com or contact your local HALO® distributor

HALO® and Fused-Core® are registered trademarks of Advanced Materials Technology, Inc.

halocolumns.com

