

Evaluation of Positively Charged Surface Phenyl-Hexyl Stationary Phase for Separation of Basic Small Molecule Drugs

William L. Miles, Barry E. Boyes, Ph.D.
Advanced Materials Technology, Inc.



HPLC 2024

July 25th, 2024

Basic Analyte Challenges in RP-LC

Achieving good peak shapes of ionized analytes by reversed-phase liquid chromatography (RP-LC) can be difficult

- Overload at low mass loads
- High asymmetry and low efficiencies
- Does not match performance of neutral or non-ionized molecules



When using low ionic strength mobile phases, like those favored in liquid chromatography-mass spectrometry (LC-MS), these effects are worsened

Neutral vs Basic Peak Shape

90Å SPP C18 2.7µm, 2.1x100mm

MP A = H₂O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

x% B, 0.50mL/min, 35C, 1.0 µL inj, 254nm

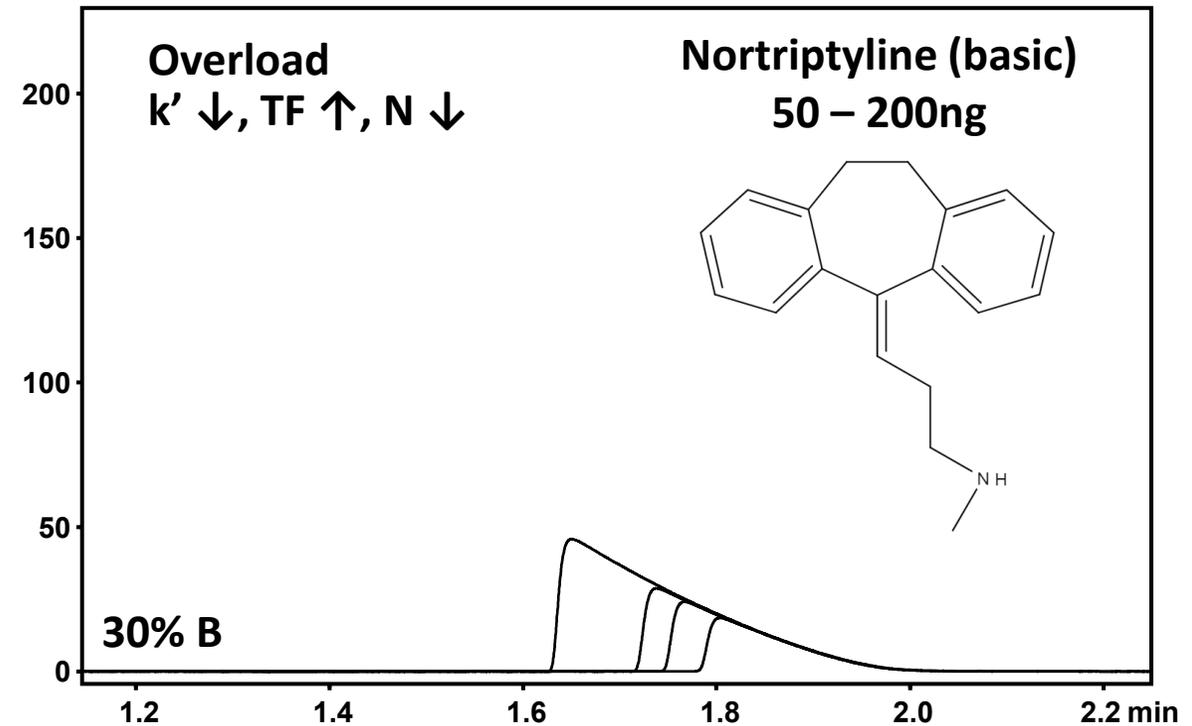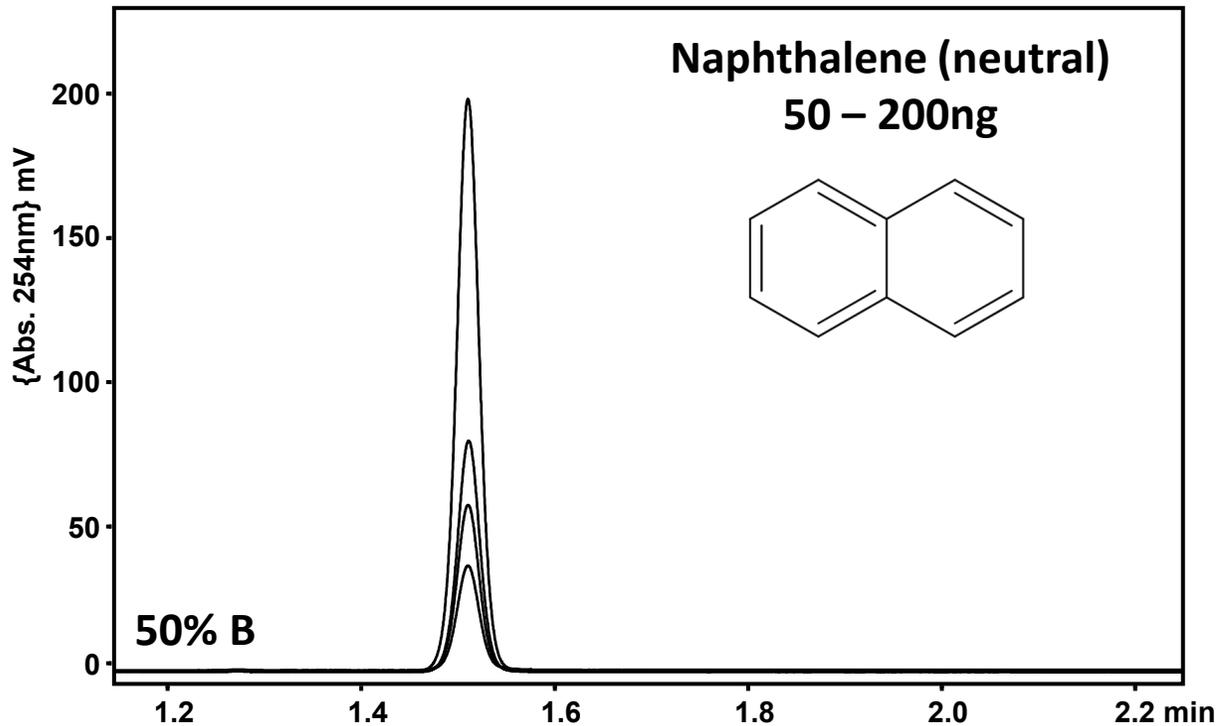

Theories on Basic Analyte Overloading

Heterogeneous adsorption of basic solutes onto stationary-phase
[1, Gritti et al. 2014]

Low Mass Loads: Adsorption onto a small number of high energy sites near the surface [1]

Higher Mass Loads: Adsorption onto a larger number of low energy sites, solute-solute repulsion [1]

Mutual repulsion of basic solutes adsorbed to bonded-phase [2, McCalley 2003]

[1] F. Gritti, G. Guiochon, Effects of the surface concentration of fixed charges in C18-bonded stationary phases on the adsorption process and on the preparative chromatography of small ionizable compounds, *Journal of Chromatography A* 1372 (2014) 42–54. <https://doi.org/10.1016/j.chroma.2014.10.003>.

[2] D.V. McCalley, Rationalization of Retention and Overloading Behavior of Basic Compounds in Reversed-Phase HPLC Using Low Ionic Strength Buffers Suitable for Mass Spectrometric Detection, *Anal. Chem.* 75 (2003) 3404–3410. <https://doi.org/10.1021/ac020715f>.

Mobile Phases Considerations

Common modifications to improve basic peak shapes in LC-UV

- Increase mobile phase pH
- Ion pairing reagents (trifluoroacetic acid)
- Increase ionic strength (ammonium formate, phosphate buffer)

LC-MS is commonly used for analysis of pharmaceuticals

- Impurities, unknowns, and metabolites

High sensitivity LC-MS (ESI) analysis require volatile, low ionic strength mobile phases for efficient ionization


- 0.1% formic acid preferred

Basic Small Molecule Pharmaceuticals


Many small molecule pharmaceuticals contain basic functional groups

Neutral and acidic pH separations (RP-LC)


- Basic groups will be protonated, positive charge

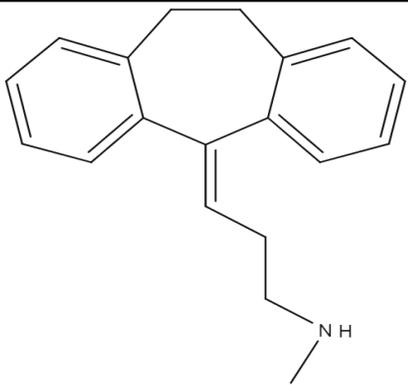

Tricyclic Antidepressants
Nortriptyline, 2°, pKa = 10.1

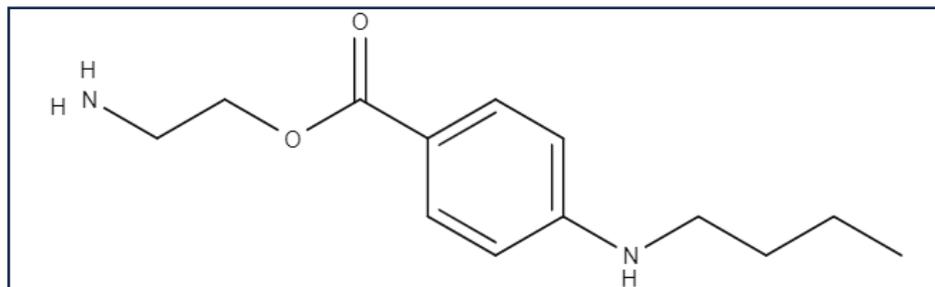

Local Anesthetics
Tetracaine, 1°, pKa = 8.5

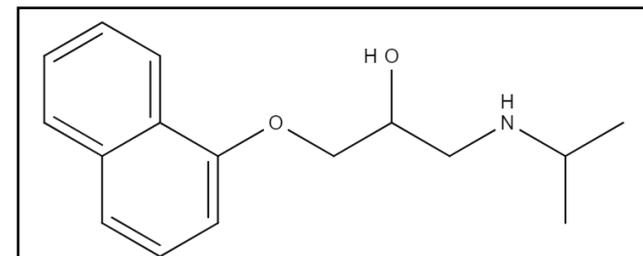

Beta-Blockers
Propranolol, 2°, pKa = 9.5

Bonded-Phase Technology

Addition of a positively charged ligand to the silica particle surface can improve basic peak shape

- Positive charge under acidic conditions
- Repulses cationic solutes, screens from negative interactions

Many versions of this technology have been described

- Commonly found with C18 bonded-phase, on FPP and SPP's


These are reversed-phase chromatographic columns

- Not ion-pairing or mixed-mode

Bonded-Phase Technology: Positively Charged Surface

90Å HALO PCS C18 2.7μm

- Superficially porous particle
- 125 m²/g surface area
- End capped

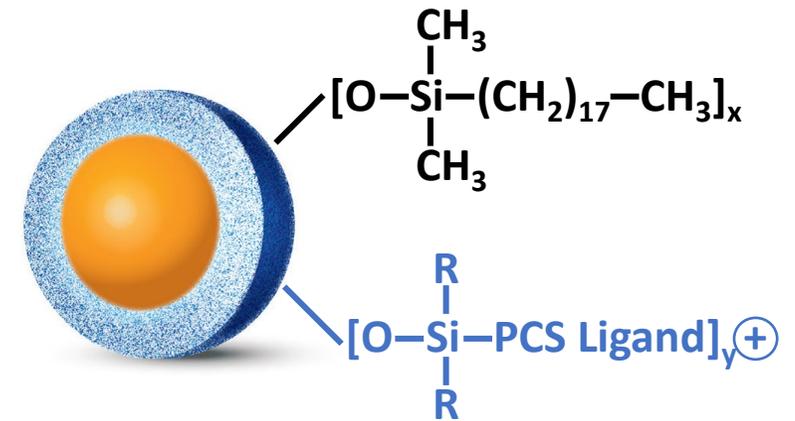

Careful selection of the PCS ligand

- Correct ligand chemistry and coverage (μmol/m²)
- [C18] > [PCS]

Designed for separations of basic small molecules, using low ionic strength mobile phases

Change in Primary Ligand: Phenyl-Hexyl

HALO[®] PCS technology: Phenyl-Hexyl

- Common RP-LC bonded-phase
- Aromatic nature, with alternative selectivity

90Å HALO PCS Phenyl-Hexyl 2.7μm



- Superficially porous particle
- 125 m²/g surface area
- End capped



Unique combination of SPP + positively charged ligand + Phenyl-Hexyl

Improvements with HALO® PCS Phenyl-Hexyl

2.1x100mm, x% B, 0.50mL/min, 35C, 1.0 µL inj, 254nm
MP A = H2O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

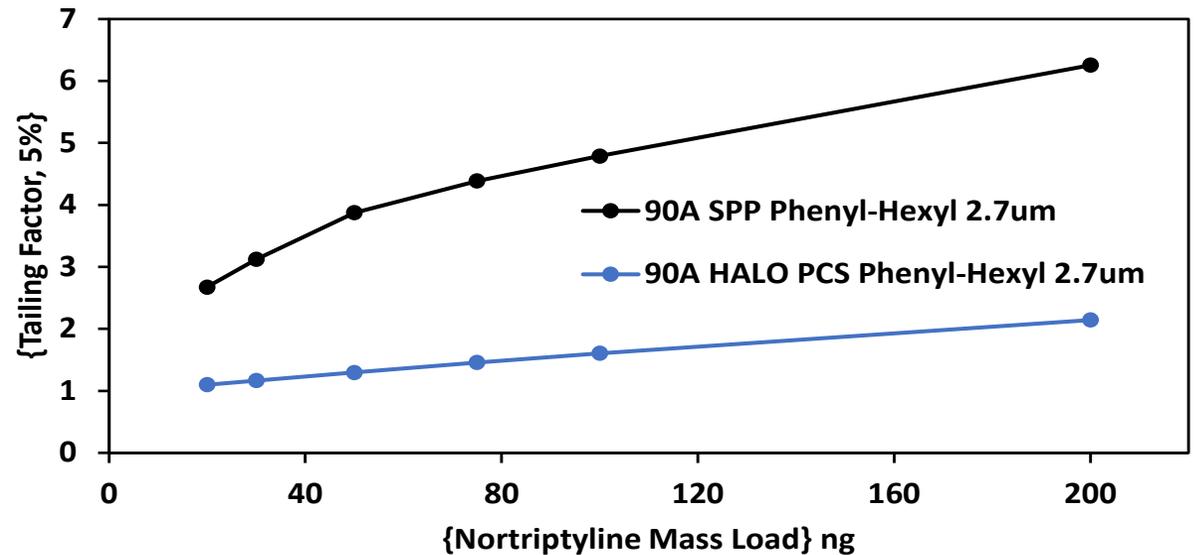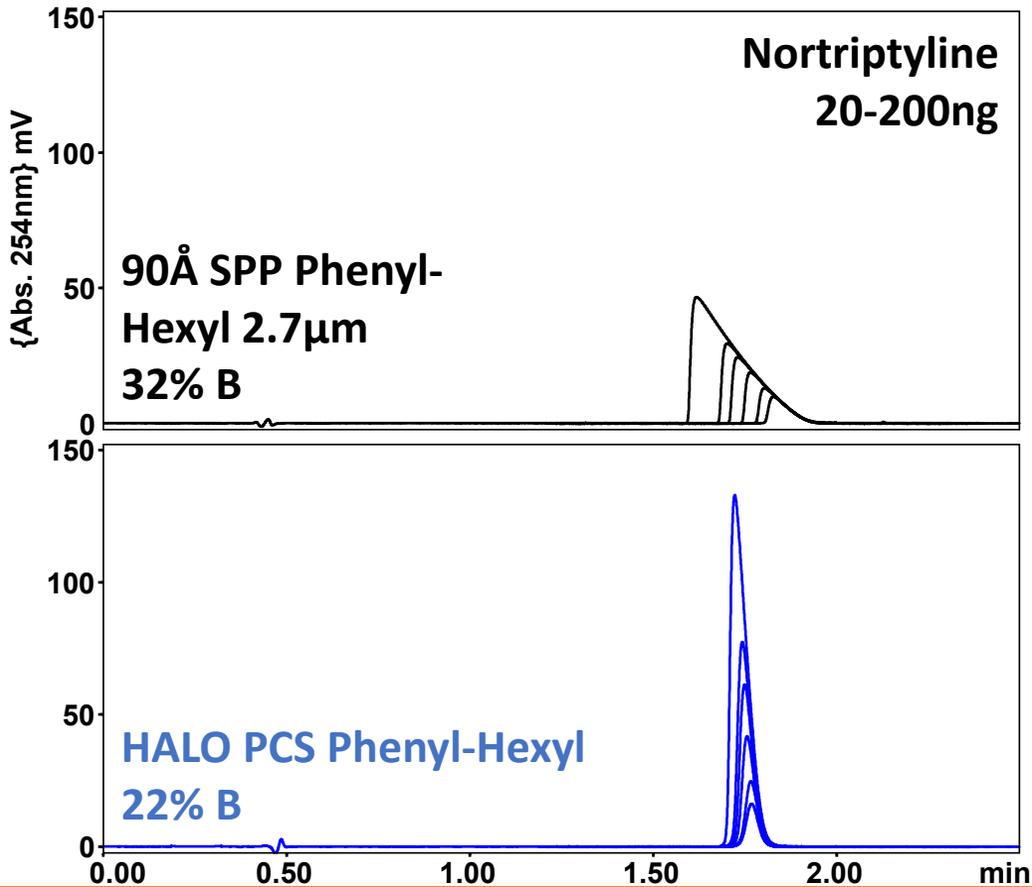

Addition of positively charged ligand:

- Improvements in peak tailing and plates
- Decrease in retention times
- PCS Phenyl-Hexyl and C18

Evaluation of HALO® PCS Phenyl-Hexyl

Separations of basic pharmaceuticals

- Differences in selectivity
- Load tolerance performance
- Imipramine applications (spiked and standard impurity)

Comparisons “In-class”


- Commercially available SPP's with C18 bonded-phase and a charged surface modification
- 90-100Å pore size, 2.7µm particle size
- 2.1x100mm columns

Mobile Phases: H₂O and ACN with 0.1% formic acid

Shimazu Nexera UHPLC PDA

Changes in Basic Analyte Selectivity


2.1x100mm, x% B, 0.50mL/min, 30C, 1.0 μ L inj, 230nm
MP A = H₂O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

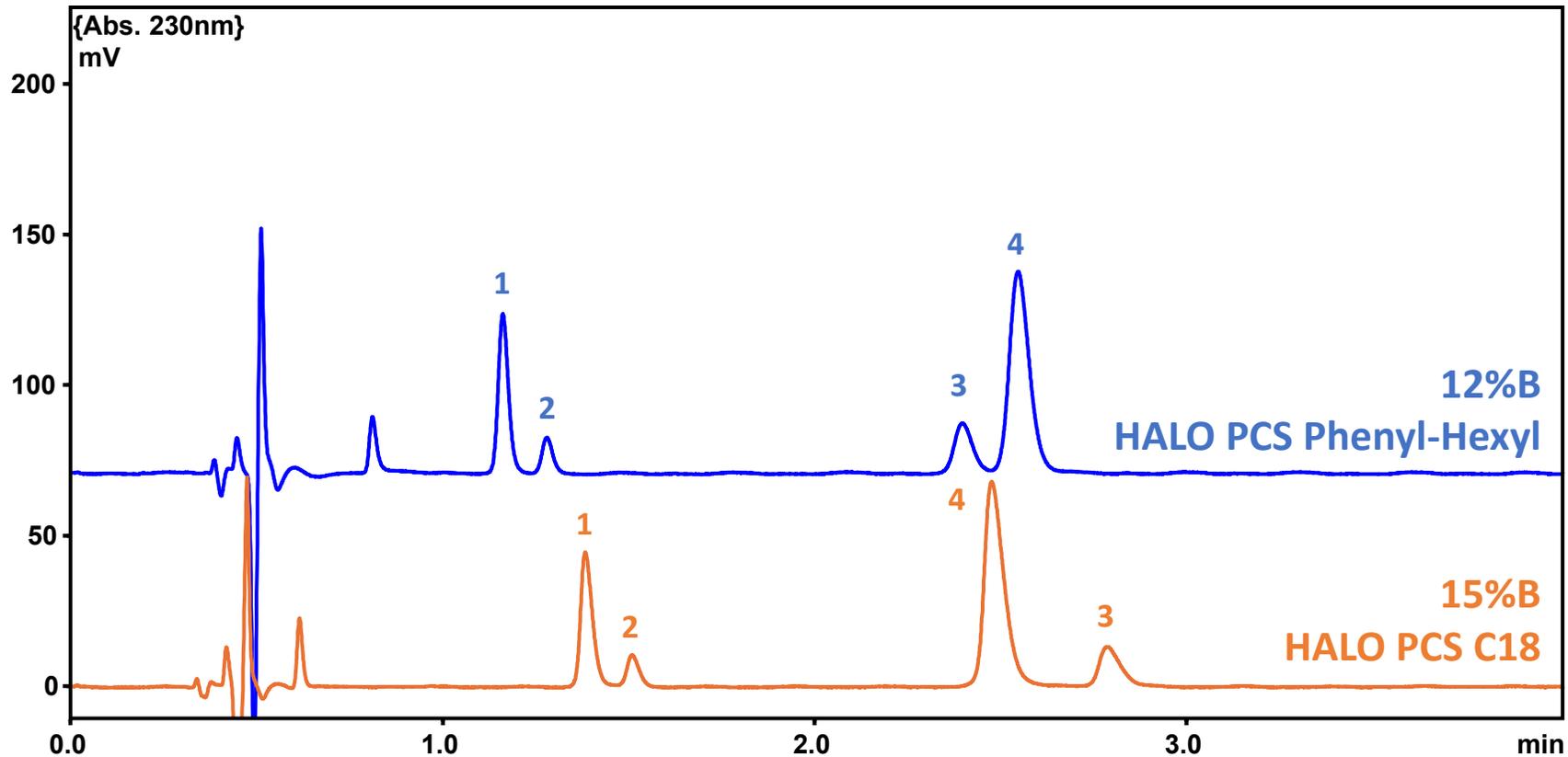


Mixture of Beta-Blockers and
Local Anesthetics
(0.030 mg/mL)

Peak Identities

1. Bisoprolol (beta-blocker)
2. Bupivacaine (local anesthetic)
3. Tetracaine (local anesthetic)

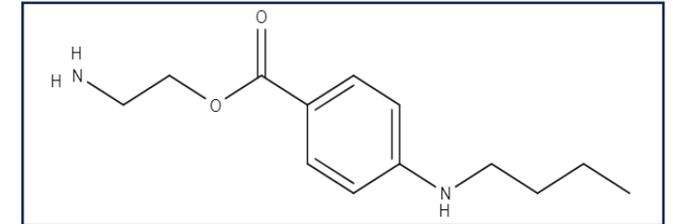


4. R-propranolol (beta-blocker)

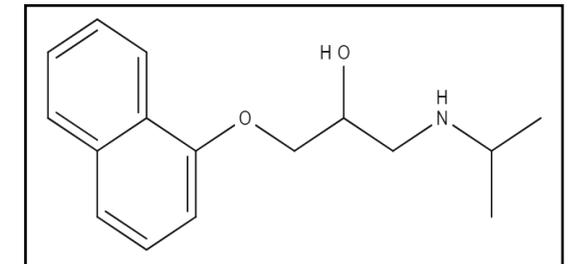

Changes in Basic Analyte Selectivity

2.1x100mm, x% B, 0.50mL/min, 30C, 1.0 μ L inj, 230nm
MP A = H₂O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

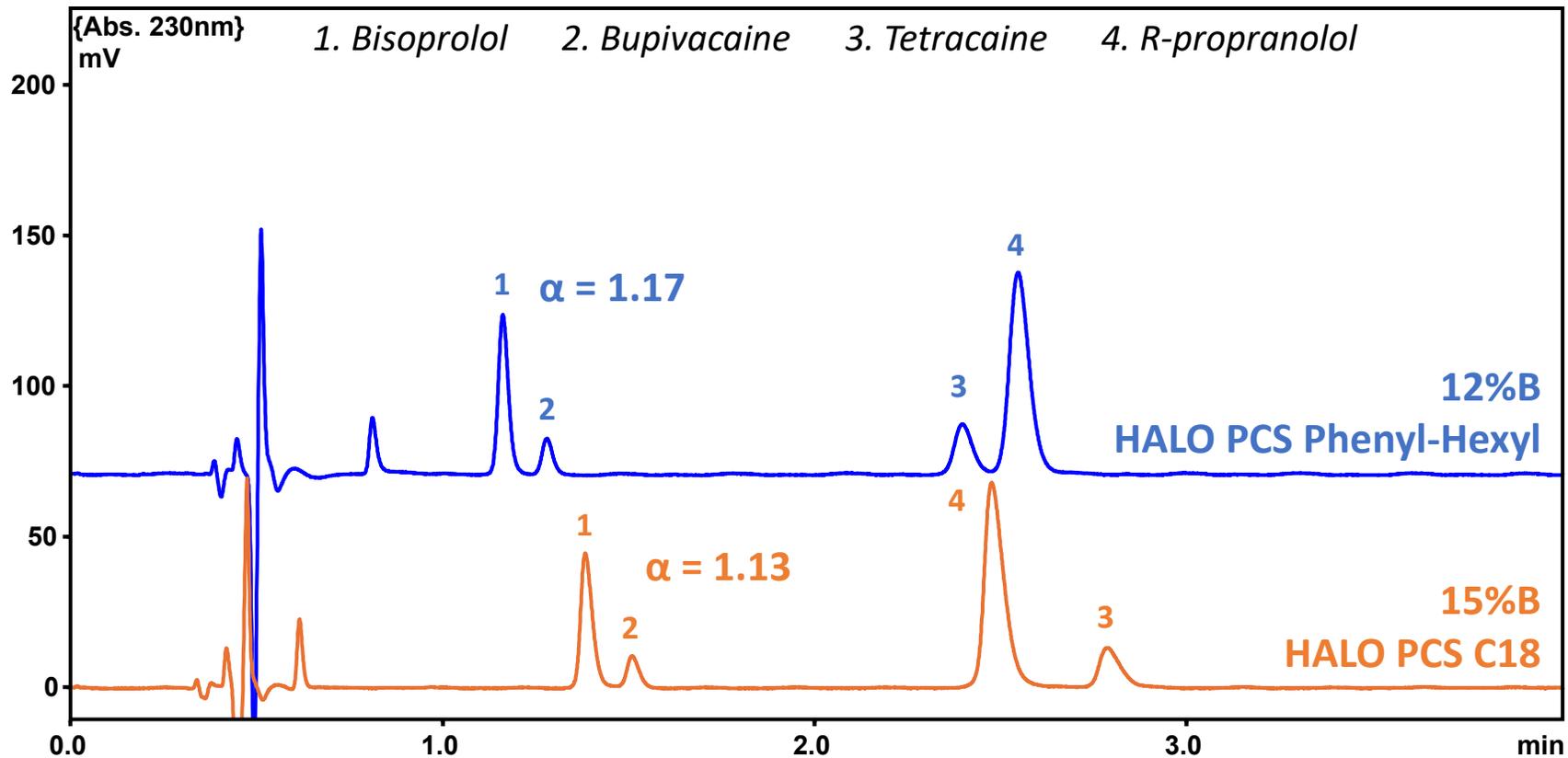

Mixture of Beta-Blockers and Local Anesthetics (0.030 mg/mL)

Retention:


- Adjusted ACN %, closer Rt match

Selectivity (α):

- Retention order switch peaks 3 & 4
- Small change for peaks 1 & 2

Changes in Basic Analyte Selectivity

2.1x100mm, x% B, 0.50mL/min, 30C, 1.0 μ L inj, 230nm
MP A = H₂O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

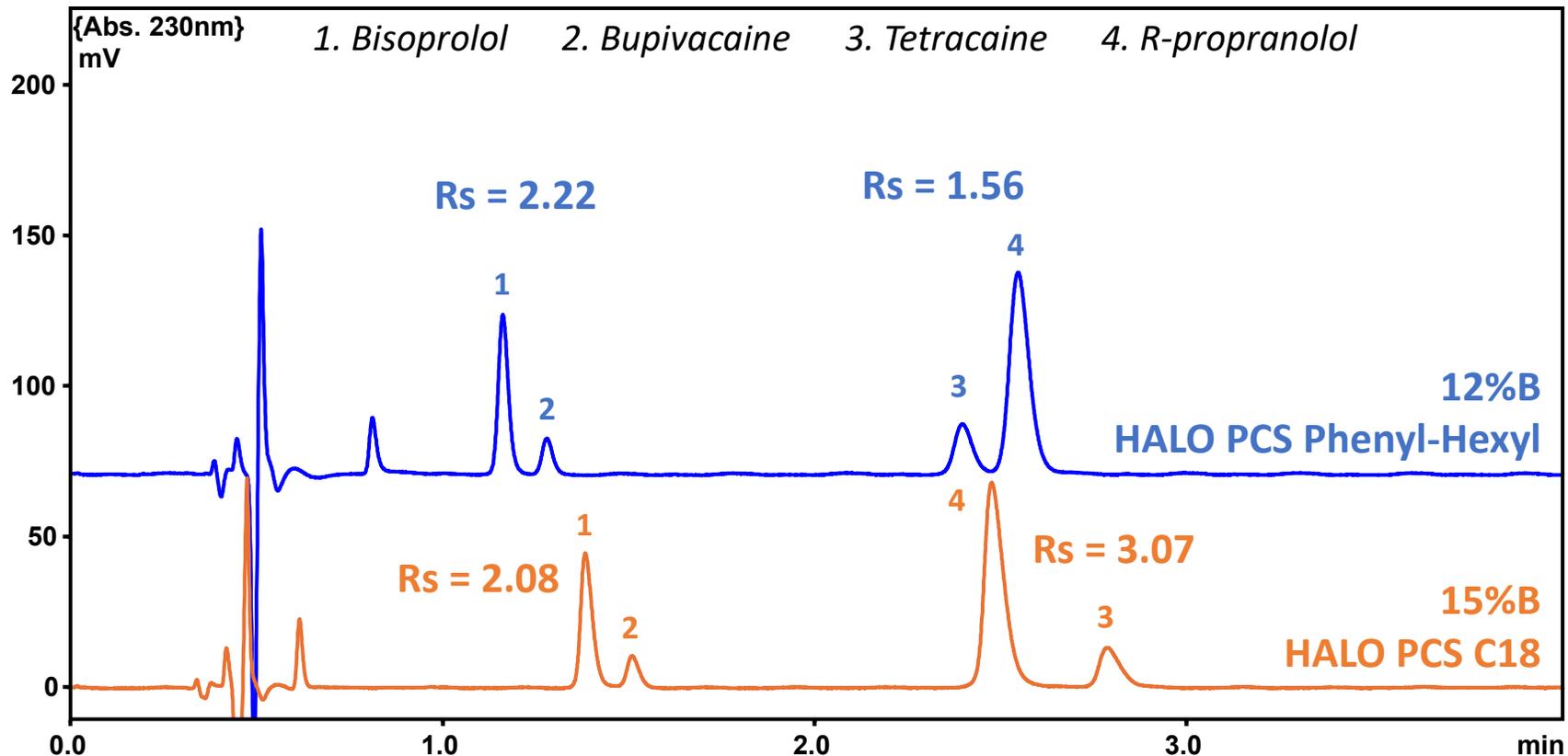


Mixture of Beta-Blockers and Local Anesthetics
(0.030 mg/mL)

Resolution (Rs):

- Baseline resolution for all pairs
- Changes in selectivity lead to changes in resolution

R-propranolol Load Tolerance

2.1x100mm, x% B, 0.50mL/min, 35C, 1.0 μ L inj, 280nm
MP A = H₂O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

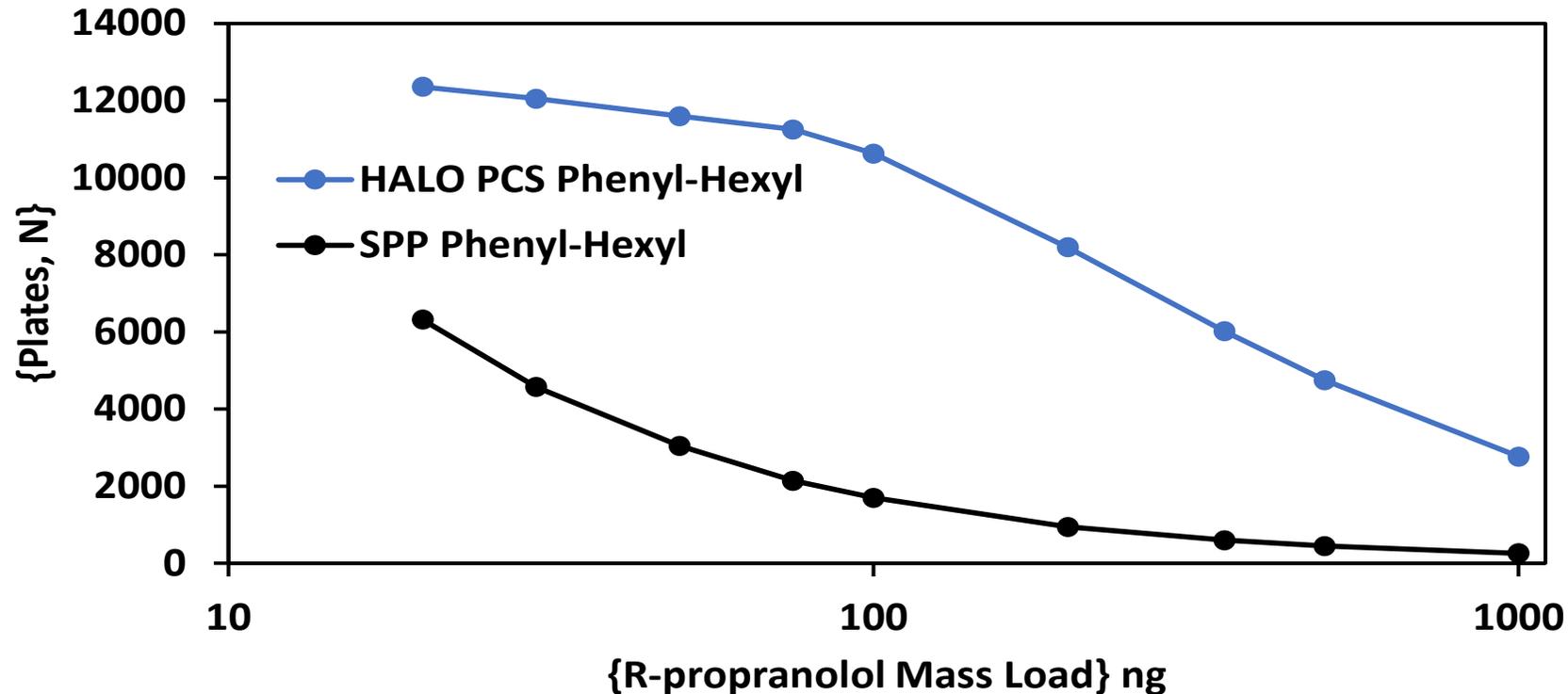

20 to 1,000 ng injected

Retention Factor, k' 20ng


- HALO[®]PCS Phenyl-Hexyl = 3.18
- SPP Phenyl-Hexyl = 3.27

7x increase in $N_{50\%}$ loss

- HALO[®]PCS Phenyl-Hexyl = 339 ng
- SPP Phenyl-Hexyl = 49 ng

R-propranolol Load Tolerance

2.1x100mm, x% B, 0.50mL/min, 35C, 1.0 μ L inj, 280nm
MP A = H₂O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

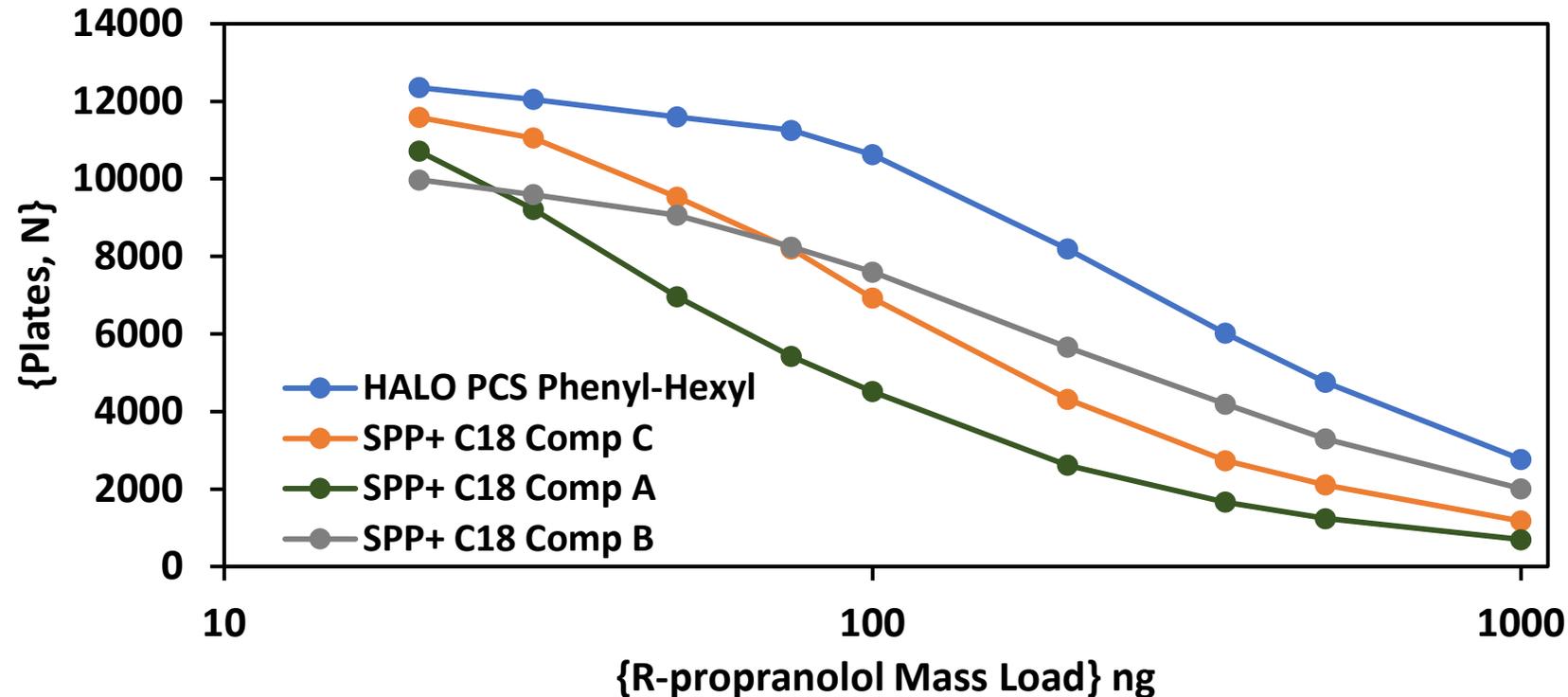

Retention factor, k' 20ng

- HALO PCS Phenyl-Hexyl = 3.18
- SPP+ C18 Comp A = 2.99
- SPP+ C18 Comp B = 3.44
- SPP+ C18 Comp C = 3.29

Overload seen for all columns as mass load increased (still ionized)

HALO PCS Phenyl-Hexyl demonstrates highest column efficiencies across entire range

Load Tolerance Comparison

2.1x100mm, x% B, 0.50mL/min, 35C, 1.0 µL inj, 280nm
MP A = H2O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

Column	R-propranolol		Nortriptyline	
	N 20ng	{N _{50%} } ng	N 20ng	{N _{50%} } ng
HALO PCS Phenyl-Hexyl	12357	339	12727	208
SPP+ C18 Comp A	10713	77	10364	63
SPP+ C18 Comp B	9972	269	9938	185
SPP+ C18 Comp C	11584	143	11868	112

Retention factors matched


Similar curves, but analyte dependence
(propranolol > nortriptyline)

HALO® PCS Phenyl-Hexyl

- Provides highest column efficiencies at all tested load
- Largest N_{50%} values for both analytes

Imipramine Spiked Impurity Analysis

2.1x100mm, 90Å HALO PCS Phenyl-Hexyl 2.7µm,
17% B, 0.50mL/min, 35C, 1.0 µL inj, 254nm
MP A = H2O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

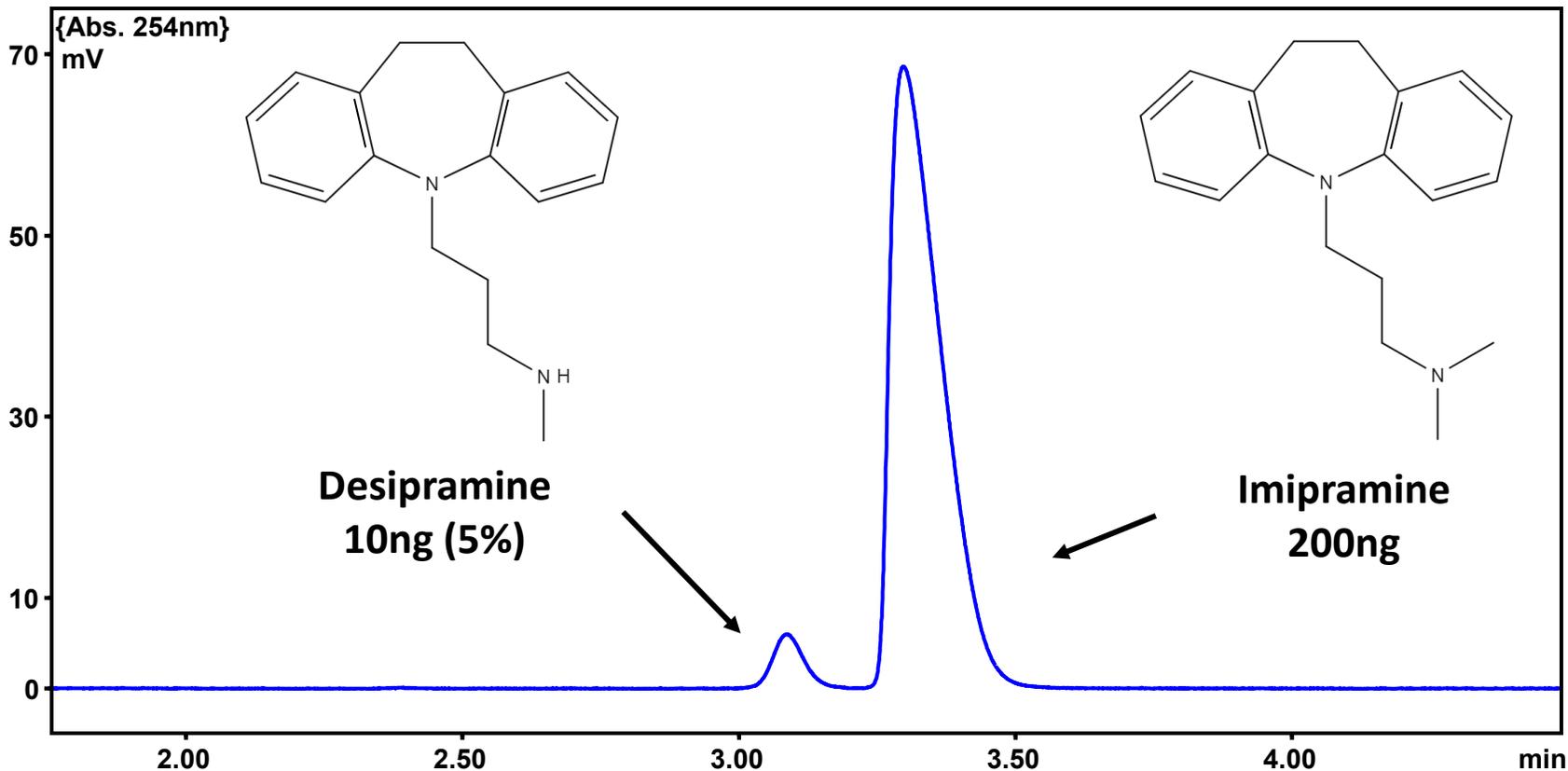

Isocratic separation of imipramine and closely related impurity


Desipramine = N-demethylation impurity

Moderately high load

- Tailing and loss of efficiency expected for imipramine

Imipramine Spiked Impurity Analysis

2.1x100mm, 90Å HALO PCS Phenyl-Hexyl 2.7µm,
17% B, 0.50mL/min, 35C, 1.0 µL inj, 254nm
MP A = H2O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

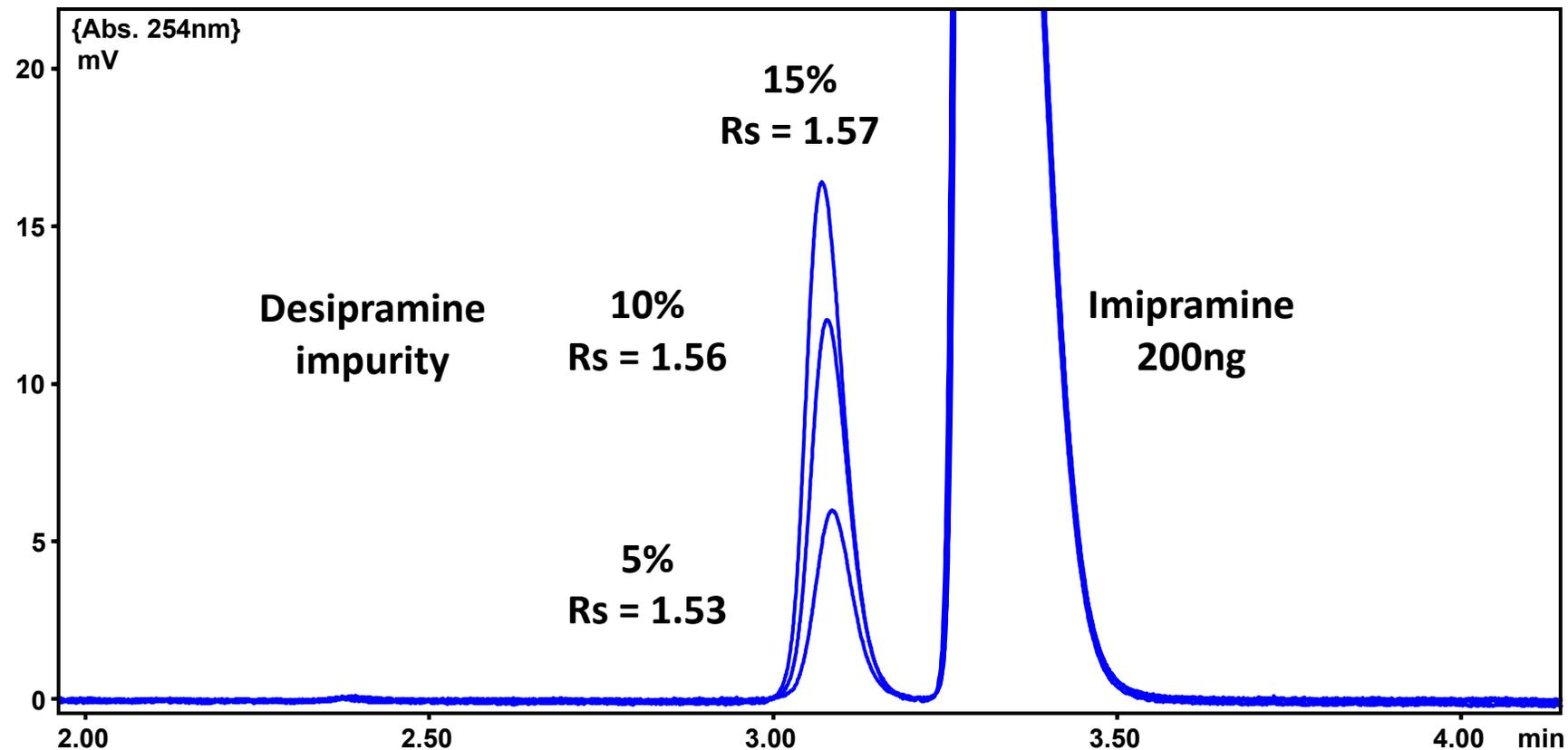


Isocratic separation (17% ACN)
of imipramine and closely
related impurity

Baseline resolution is
maintained as impurity
concentration increased

Advantage of improved column
efficiencies over a range of
mass

Imipramine Impurity Analysis LC-PDA

2.1x100mm, 90Å HALO PCS Phenyl-Hexyl 2.7µm,
16% B, 0.50mL/min, 35C, 2.0 µL inj, 254nm
MP A = H2O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

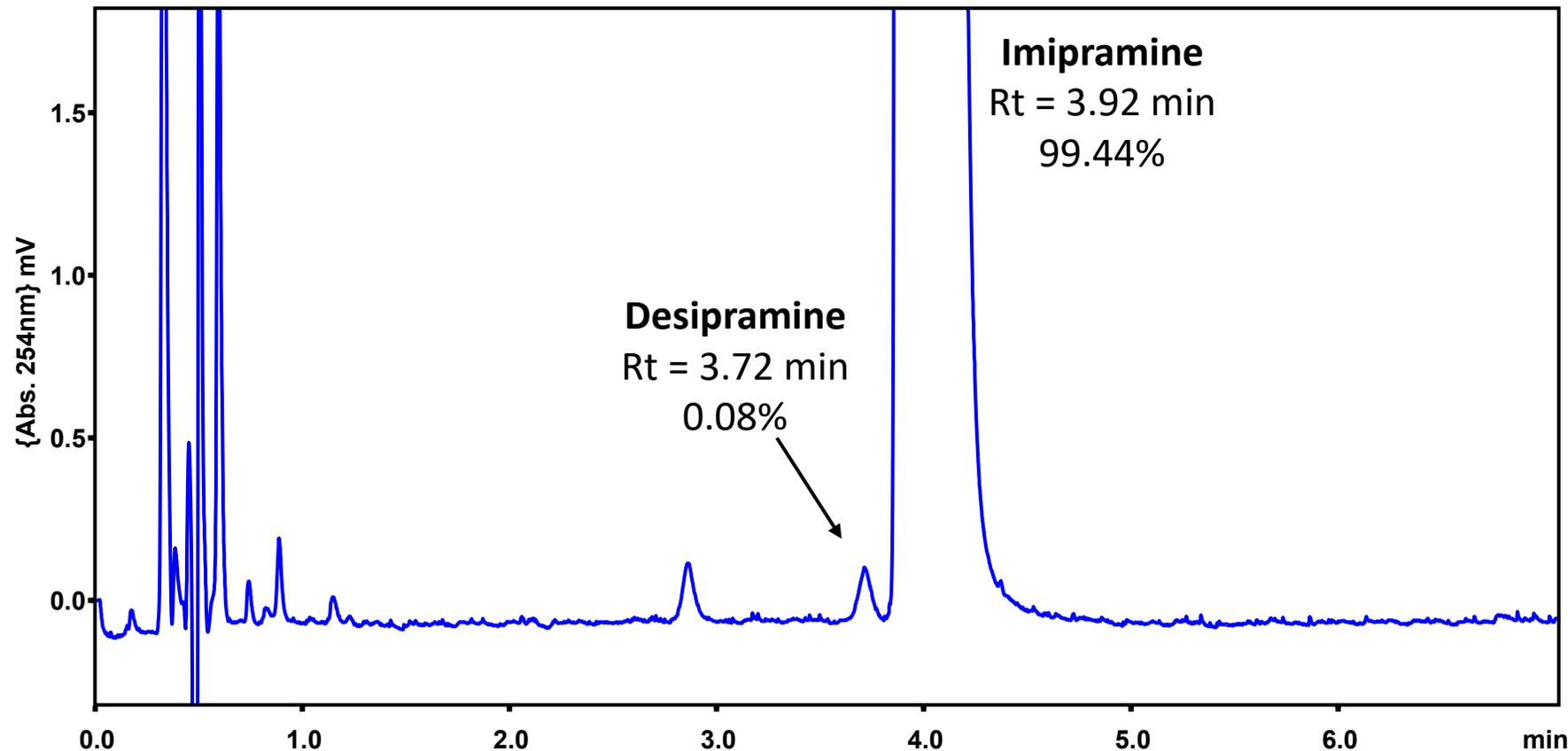

LC-PDA Isocratic Separation

400ng inj. imipramine standard

Percent Area


- 99.44% imipramine
- 0.56% impurities (0.01 – 0.33 %)
- Desipramine identified

Improved load tolerance allows for detection of trace impurities

Imipramine Impurity Analysis LC-MS

2.1x100mm, 90Å HALO PCS Phenyl-Hexyl 2.7µm,
10-25% B in 10min, 0.40mL/min, 35C, 2.0 µL inj,
MP A = H2O + 0.1% formic acid, MP B = ACN + 0.1% formic acid

Thermo Q Exactive Orbitrap HF, ESI: 3.3 kV, Tcap 320 C, Rs 45,000,
50 ms; ddMS² NCE 30%, Rs 30,000, 50 ms

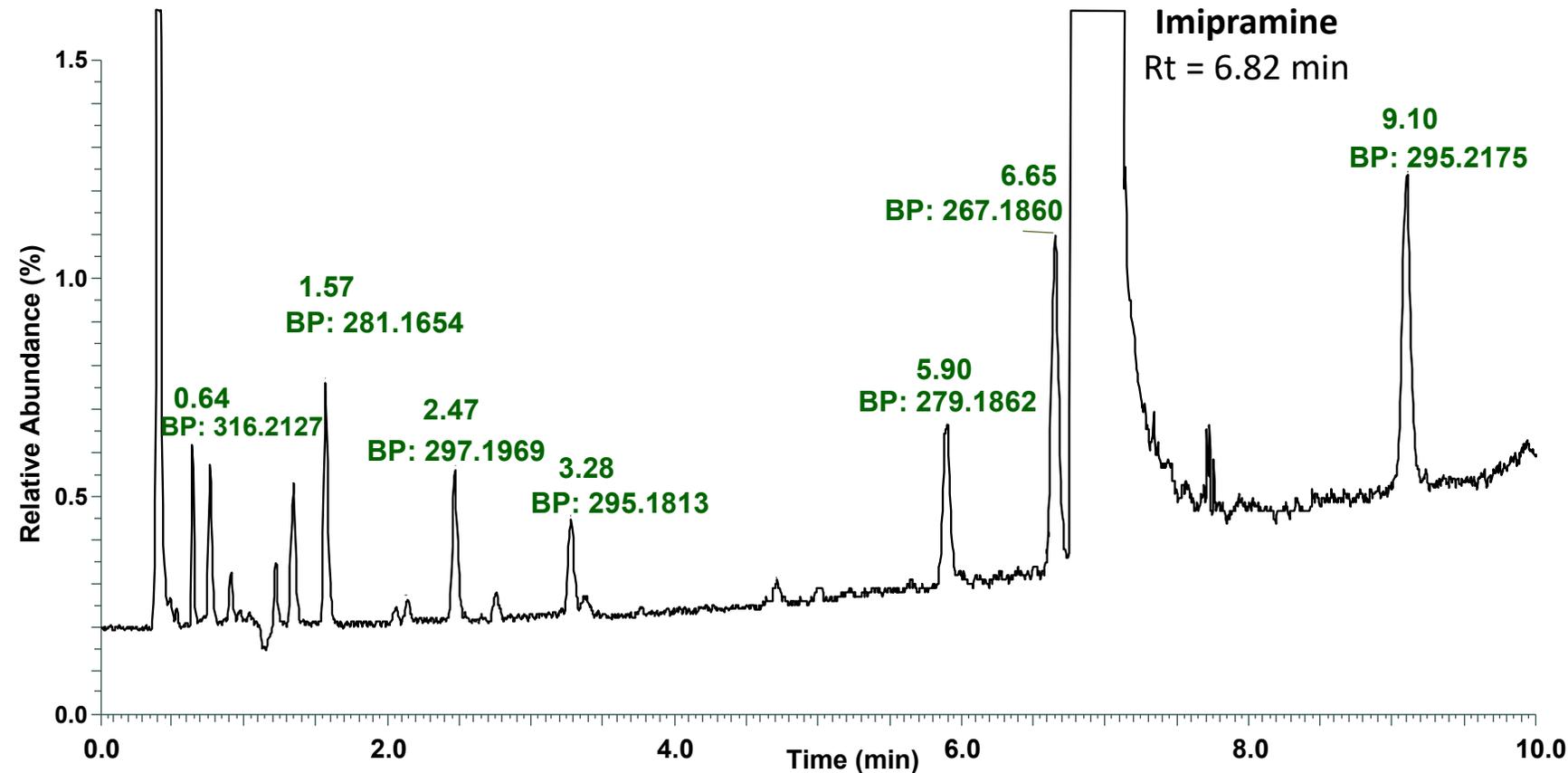

LC-MS Gradient elution method
(400ng)

Impurities \leq 1% rel. abundance

High resolution accurate mass


- Measuring exact mass
- Formula match impurities

MS² Experiment

- Structural details of impurities


Imipramine Impurity Analysis LC-MS

Peak	{Rt} min	Measured m/z	Formula Match	Functional Group	MS ² Confirmation
Imipramine	6.86	281.2009	C ₁₉ H ₂₅ N ₂	[Imip]	Y
1-4	0.76 – 1.57			No TCA	N
5	2.47	297.1969	C ₁₉ H ₂₅ N ₂ O	[Imip] +O, -OH alkyl	Y
6	3.27	295.1815	C ₁₉ H ₂₃ N ₂ O ₁	No TCA	N
7	5.90	279.1863	C ₁₉ H ₂₃ N ₂	[Imip] -2H, C=C alkyl	Y
Desipramine	6.65	267.1860	C ₁₈ H ₂₃ N ₂	[Imip] – CH ₃ , N-methyl	Y
8	9.10	295.2175	C ₂₀ H ₂₇ N ₂	[Imip] + CH ₂ (butylamine)	Y

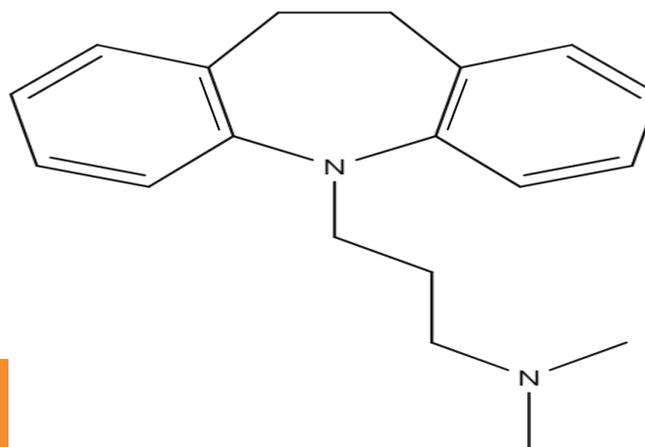

Imipramine Impurity Analysis LC-MS

Peak	{Rt} min	Measured m/z	Formula Match	Functional Group	MS ² Confirmation
Imipramine	6.86	281.2009	C ₁₉ H ₂₅ N ₂	[Imip]	Y
1-4	0.76 – 1.57			No TCA	N
5	2.47	297.1969	C ₁₉ H ₂₅ N ₂ O	[Imip] +O, -OH alkyl	Y
6	3.27	295.1815	C ₁₉ H ₂₃ N ₂ O ₁	No TCA	N
7	5.90	279.1863	C ₁₉ H ₂₃ N ₂	[Imip] -2H, C=C alkyl	Y
Desipramine	6.65	267.1860	C ₁₈ H ₂₃ N ₂	[Imip] – CH ₃ , N-methyl	Y
8	9.10	295.2175	C ₂₀ H ₂₇ N ₂	[Imip] + CH ₂ (butylamine)	Y

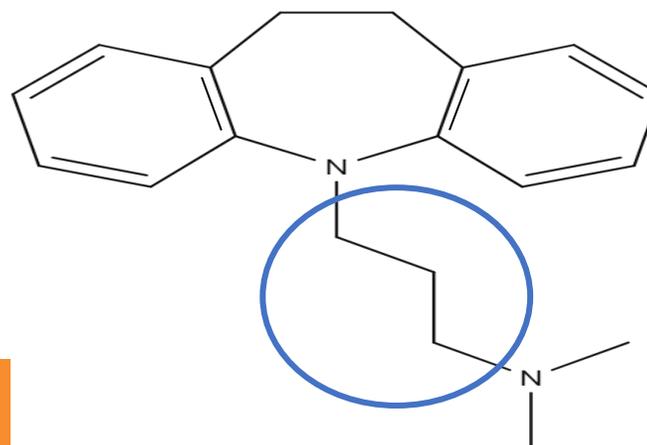

Conclusions

Changes in selectivity for HALO[®] PCS Phenyl-Hexyl compared to in-class C18's

- Another versatile option for separations

Improvements in column efficiencies over a large mass range for basic small molecule pharmaceuticals

- Phenyl-Hexyl + positively charged surface appears to be favorable for separations of these aromatic, basic pharmaceuticals
- Impurities resolution and trace impurities

Addition of positively charged ligand to the silica surface greatly improves basic peak shape

- LC-MS preferred mobile phases

Acknowledgements

Research and Development Team

- Dr. Joseph DeStefano
- Timothy Langlois
- Mark Haynes

Applications


- Dr. Stephanie Schuster
- Conner McHale

Engineering Team

- Aggie Cesbron
- Kristen O'Donnell
- Matt Jackson
- Bryant McKnight

Thank you for your time today!

Questions

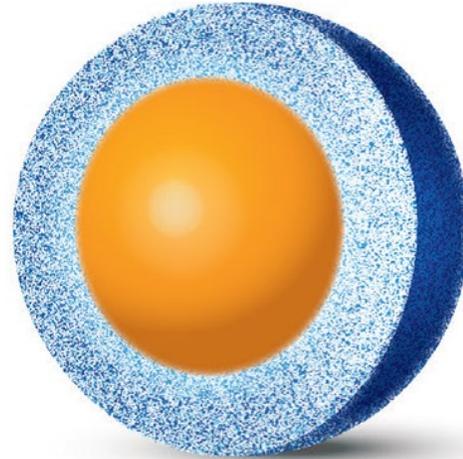



advancedmaterialstechnology



halocolumns.com



3521 Silverside Road, Suite 1-K
Quillen Building
Wilmington, DE 19810



(302) 992-8060

ISO 9001 QMS certified
All operations in Wilmington, DE
Global distribution

