

Assessment of Stationary Phase Ligands on Positively Charged Particles by an Anionic Exchange Parameter

William Miles, Stephanie Schuster, Barry Boyes, Joseph DeStefano

Advanced Materials Technology, Inc.

HPLC 2026
June 8th




advancedmaterialstechnology

Cation Modified Stationary Phases

Further our understanding of the properties of cation-modified reversed-phase (RP) column packing materials to develop high performance LC and LC-MS applications.


- Combination of primary RP ligand with a low quantity of a positively charged surface modification
 - Many commercially available options
- Improve ionized basic solute peak shape and performance, while operating in 0.1 % (v/v) formic acid mobile phases
 - High sensitivity LC-MS applications

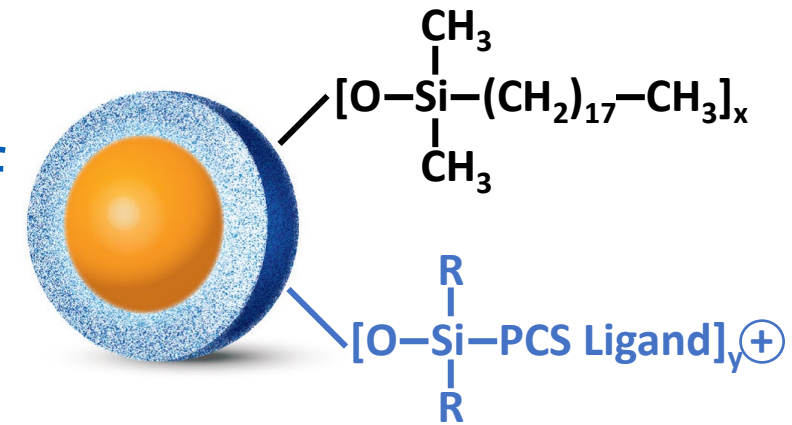

90Å HALO® PCS C18
2.7µm


Ionized Basic Solute Overload

4.6x100 mm, 1.20 mL/min, 30° C, 4.0 µL Inj, 254 nm
MP - ACN/H₂O + 0.1 % formic acid (v/v),
Amitriptyline 3 to 100 ng

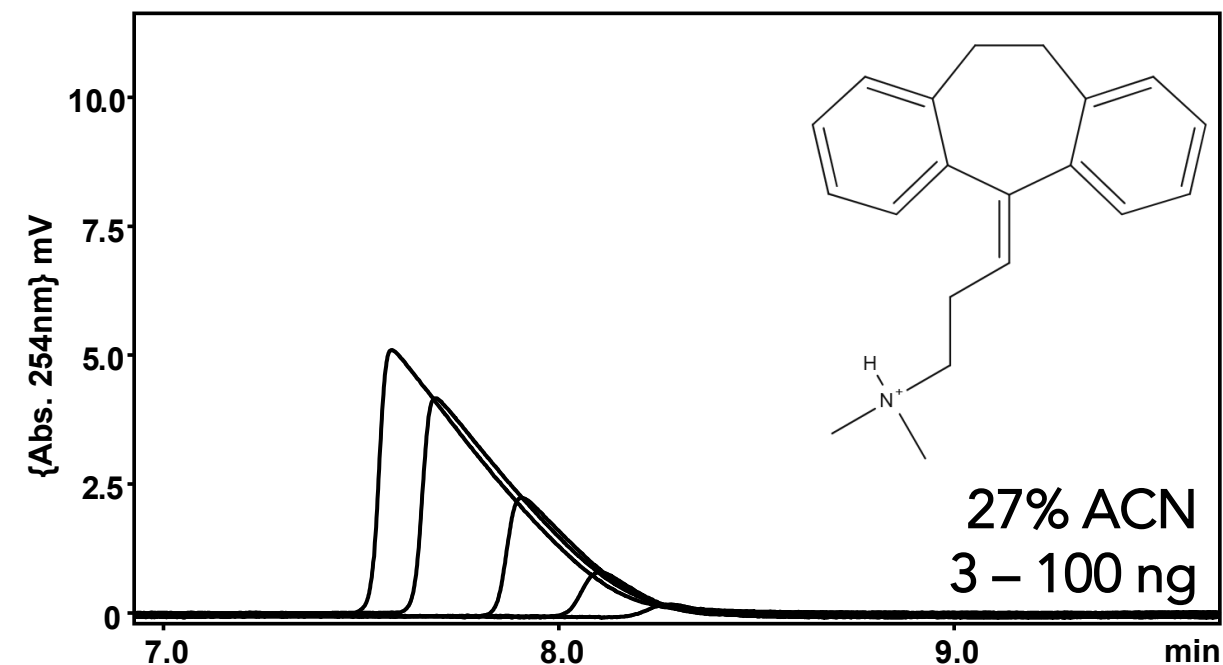


"Neutral" - 90Å C18 2.7µm SPP

Greater resistance to
ionized basic solute overload

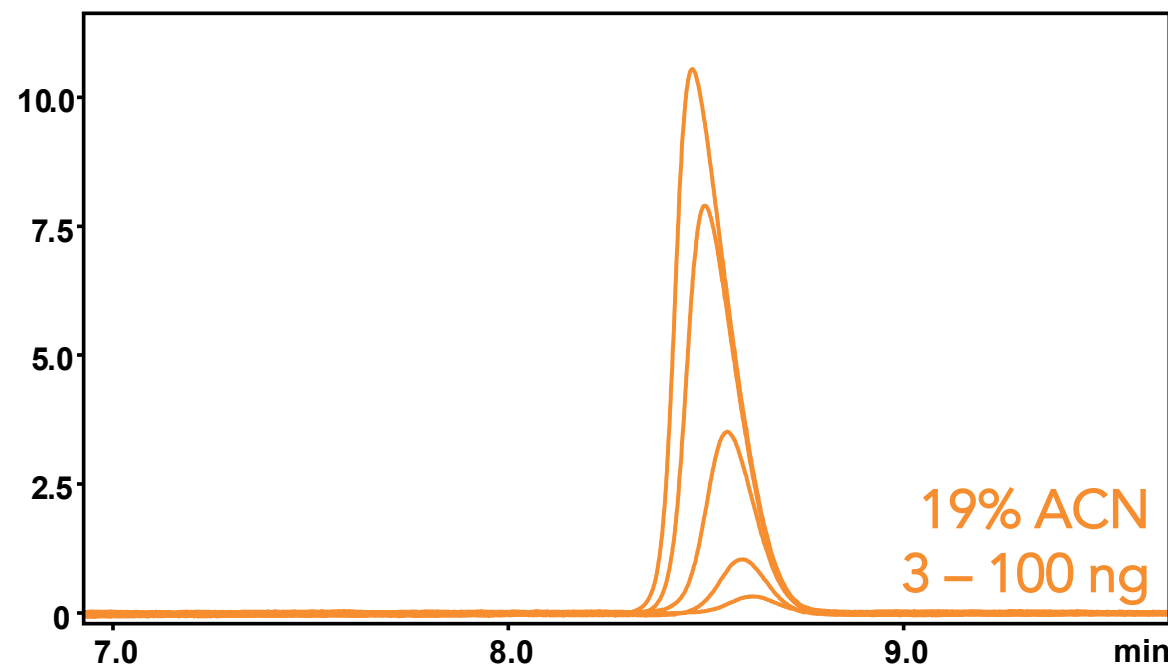

Positively Charged - 90Å PCS C18 2.7µm

Ionized Basic Solute Overload

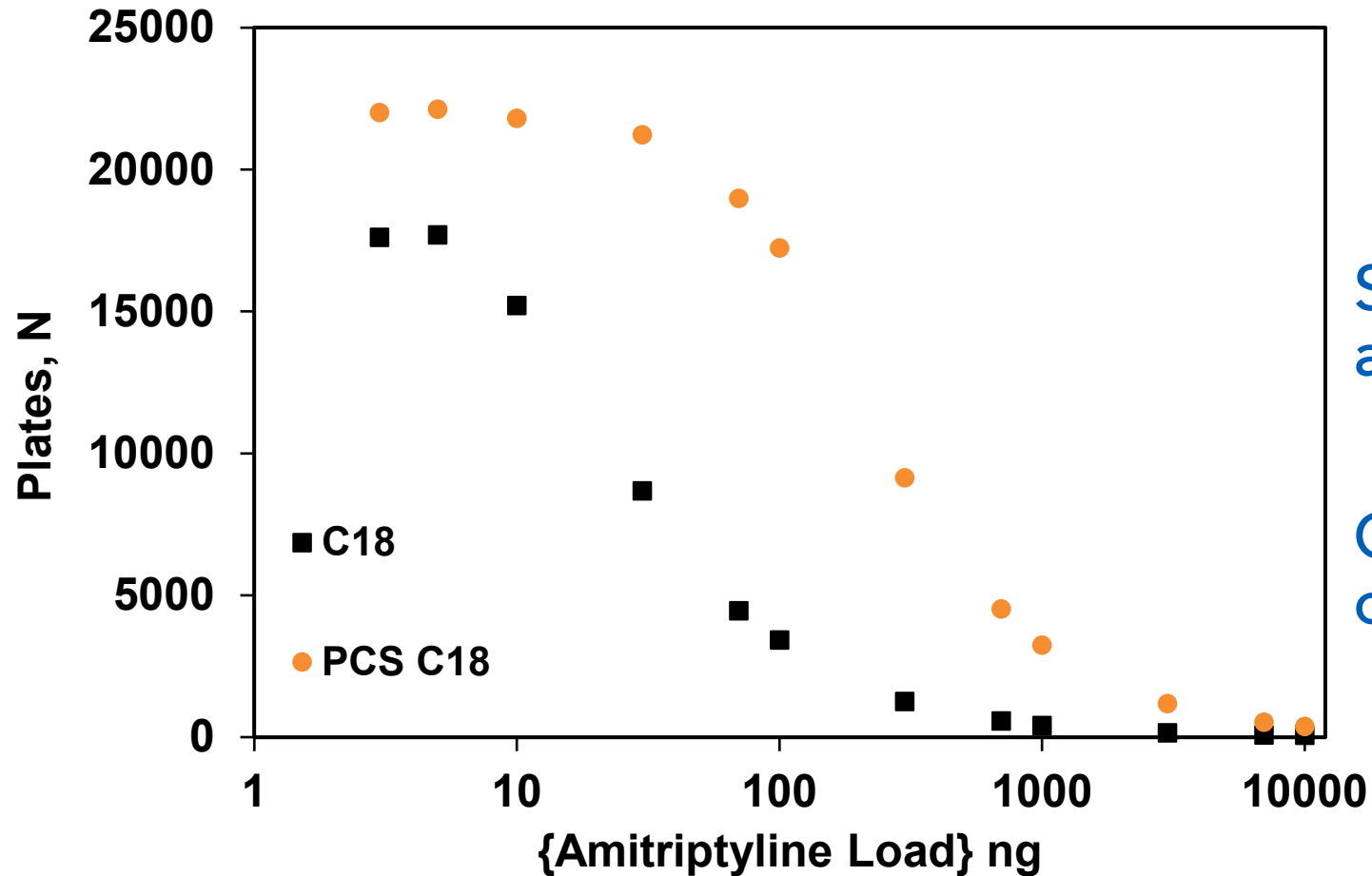


Similar trends for k'
and TF

Overload still
occurring

4.6x100 mm, 1.20 mL/min, 30° C, 4.0 μ L Inj, 254 nm, 27 and 19 % ACN, MP - ACN/H₂O + 0.1 % F.A. (v/v)

Effect of Primary Ligand

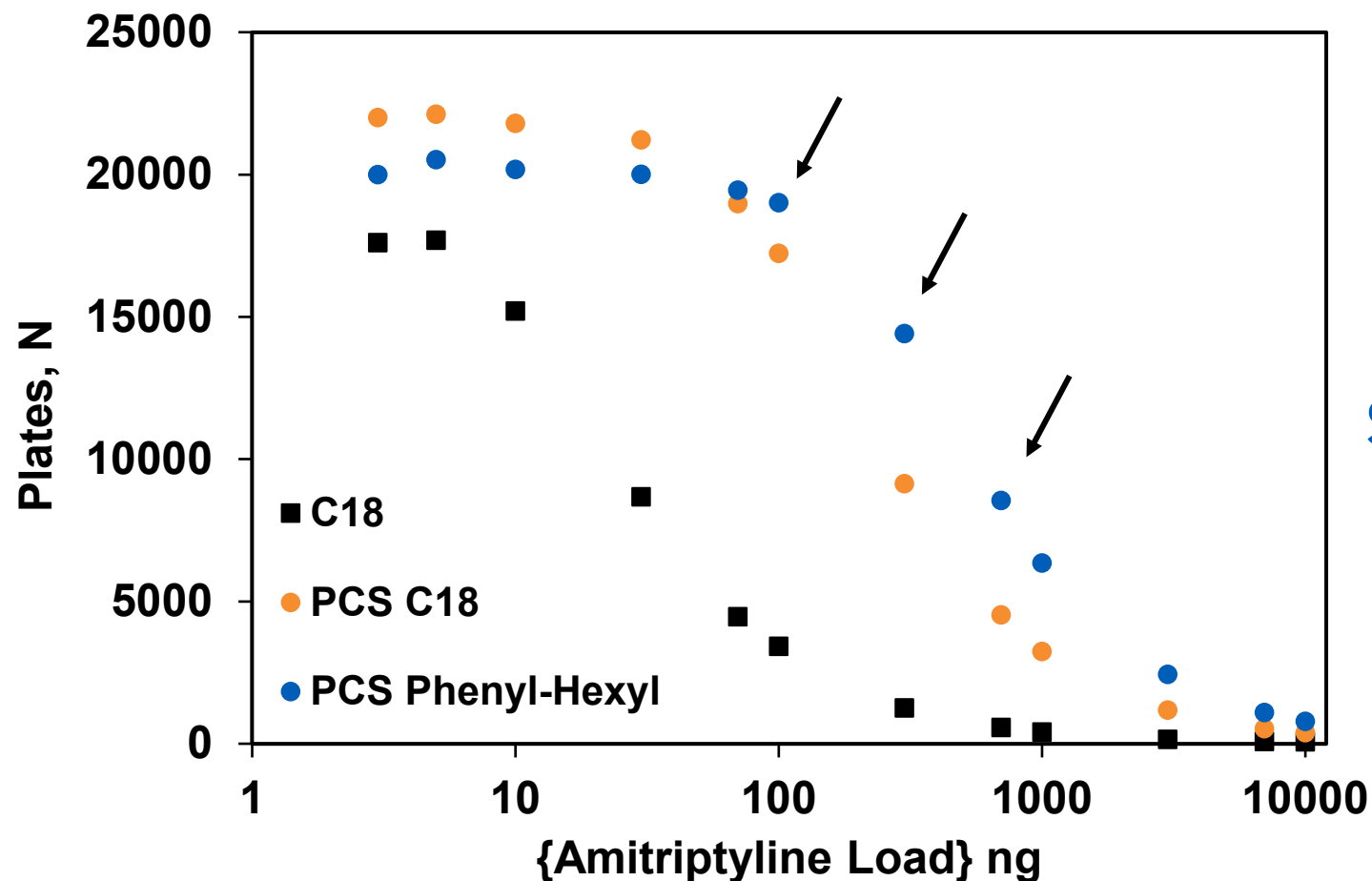

PCS Phenyl-Hexyl -
Similar improvements
in k' and TF

4.6x100 mm, 1.20 mL/min, 30°C, 4.0 μ L Inj, 254 nm, 27 / 19 / 16 % ACN, MP - ACN/H₂O + 0.1 % F.A. (v/v)

Anionic Exchange Parameter

A 2024 publication by Jolliffe et. al. described an updated anionic exchange (AE) parameter, to better discriminate between the positive character of neutral and cation modified stationary phases [1].

- Tanaka Plots – New AE Parameter $\alpha_{4\text{-OBSA} / \text{Tol}}$ [1]
 - Replace α benzenesulfonic acid (BSA) with α 4-n-octylbenzenesulfonic acid (4-OBSA)
 - Measure selectivity of anion with toluene (Tol)
- Classified the positive character of phases into three groups: neutral, low to medium, and significant [1]

[1] S. Jolliffe, J.K. Field, M.R. Euerby, P. Petersson, Characterization of reversed-phase liquid chromatographic columns containing positively charged functionality, *Journal of Chromatography A* 1730 (2024) 465059.

Anionic Exchange Parameter

What can we learn from this new AE Parameter $\alpha_{4\text{-OBSA} / \text{Tol}}$ to help in developing stationary phases?

Measure $\alpha_{4\text{-OBSA} / \text{Tol}}$ for cation modified stationary phases bonded with different primary hydrophobic ligands

- Does the primary hydrophobic ligand play a role in determining the positive character of these phases?


Design of Experiments

90Å HALO® PCS 2.7µm (SPP)


- Silica, 125 m²/g surface area
- C18 and Phenyl-Hexyl commercially available
- Controlled experiment
 - Surface area and coverage of PCS

Experiments


- $\alpha_{4\text{-OBSA/Tol}}$ AE measurements
- Ionized solute overload studies
- LC-MS Beta Blocker Separation

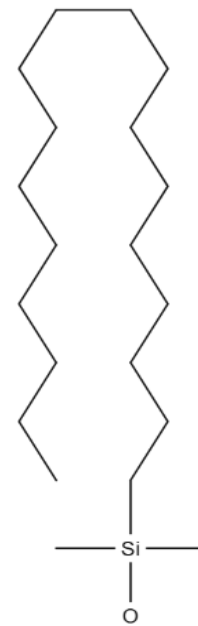


C18

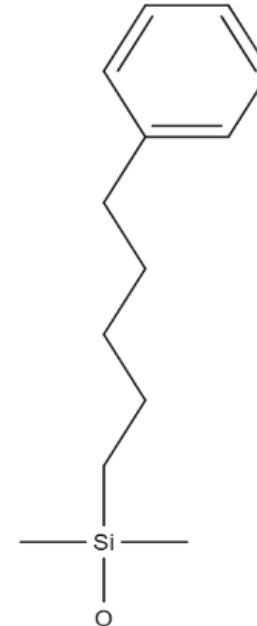


Phenyl-Hexyl

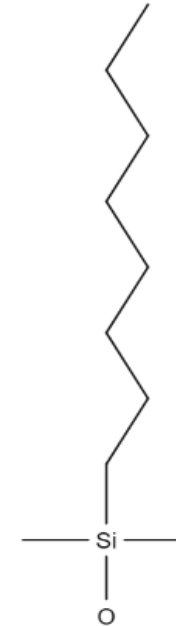

C8

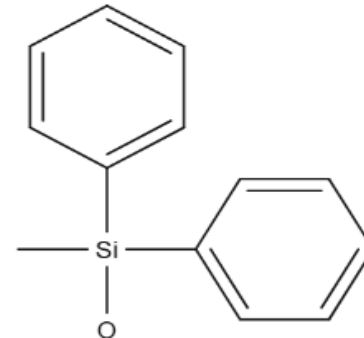


Diphenyl

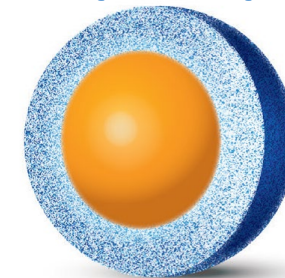

Anionic Exchange Measurement

4.6x100 mm, 1.80 mL/min, 40° C, 1.0 µL Inj, 215 nm
MP – 50/50 ACN/H₂O + 5mM KH₂PO₄ pH = 2.5 , Ref [1]
0.50 mg/mL

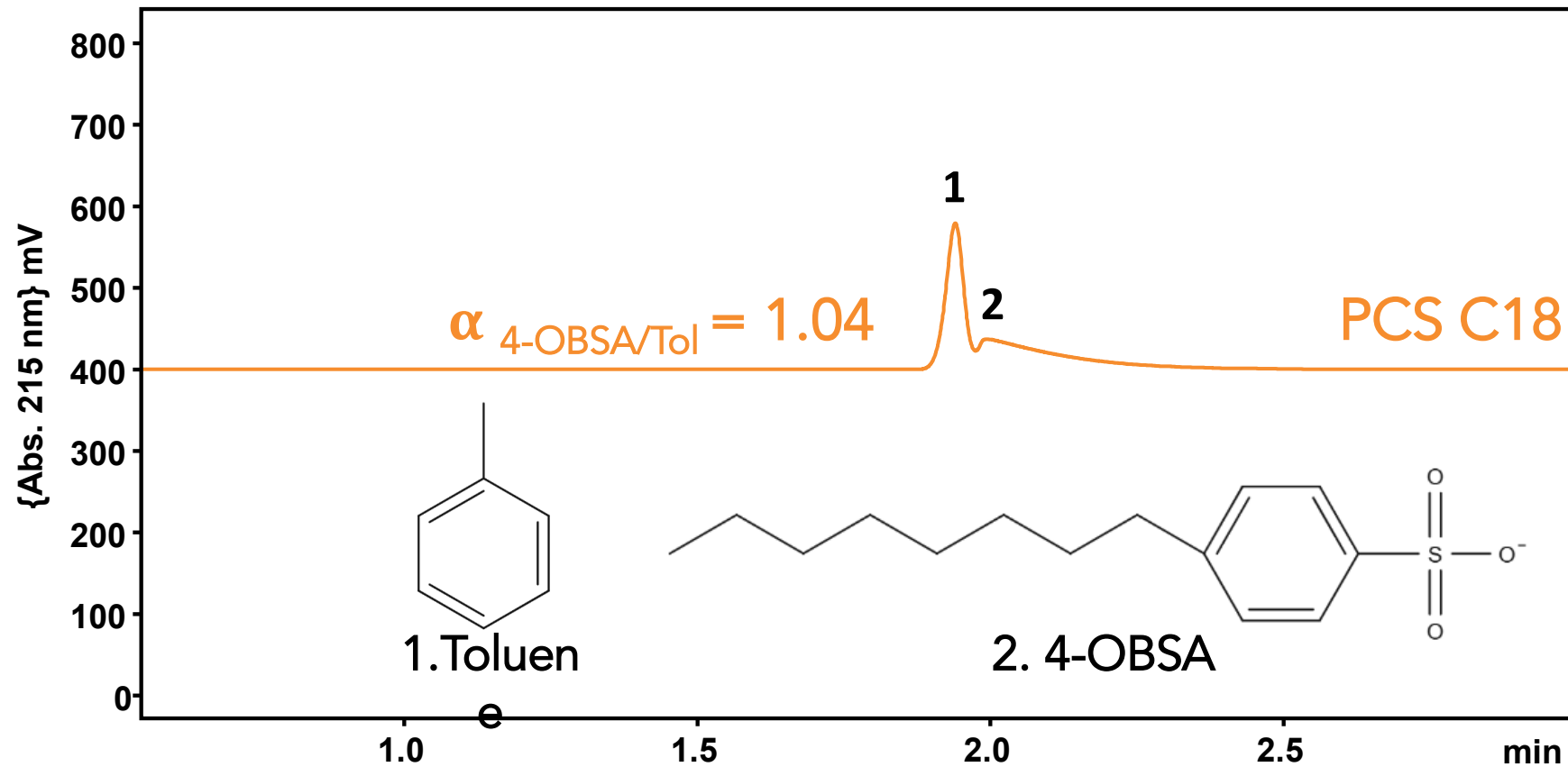


Anionic Exchange Measurement

4.6x100 mm, 1.80 mL/min, 40° C, 1.0 μ L Inj, 215 nm
MP – 50/50 ACN/H₂O + 5mM KH₂PO₄ pH = 2.5 , Ref [1]

1.Toluene (0.50 mg/mL)

2. 4-OBSA (0.50 mg/mL)

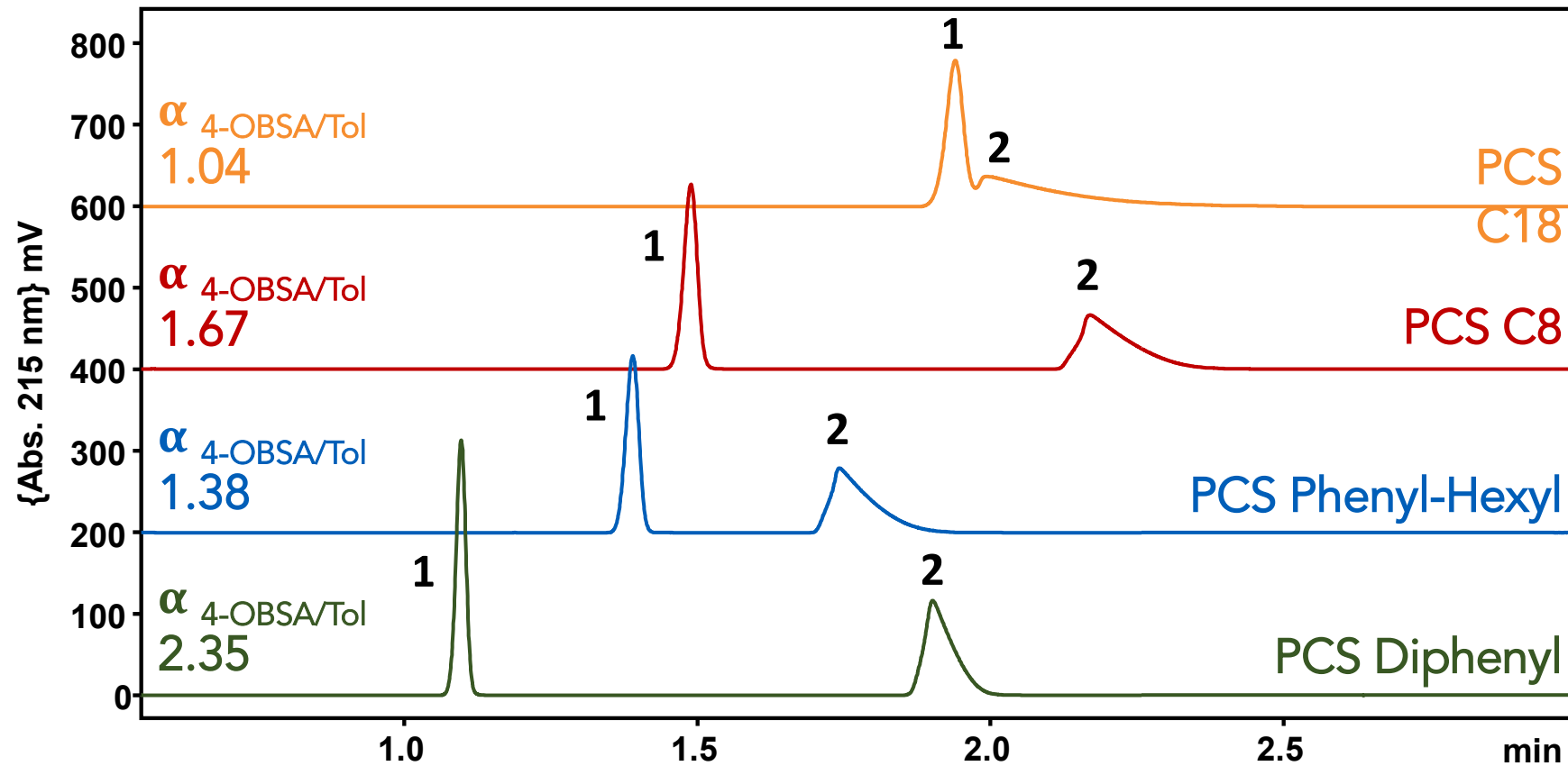


Anionic Exchange Formic Acid

Formic acid mobile phase

- Decrease in ionic strength
- 3-to-4-fold increase in α_{4-} OBSA/Tol

Positive Character of PCS Phases


Diphenyl > C8 > Phenyl-Hexyl > C18

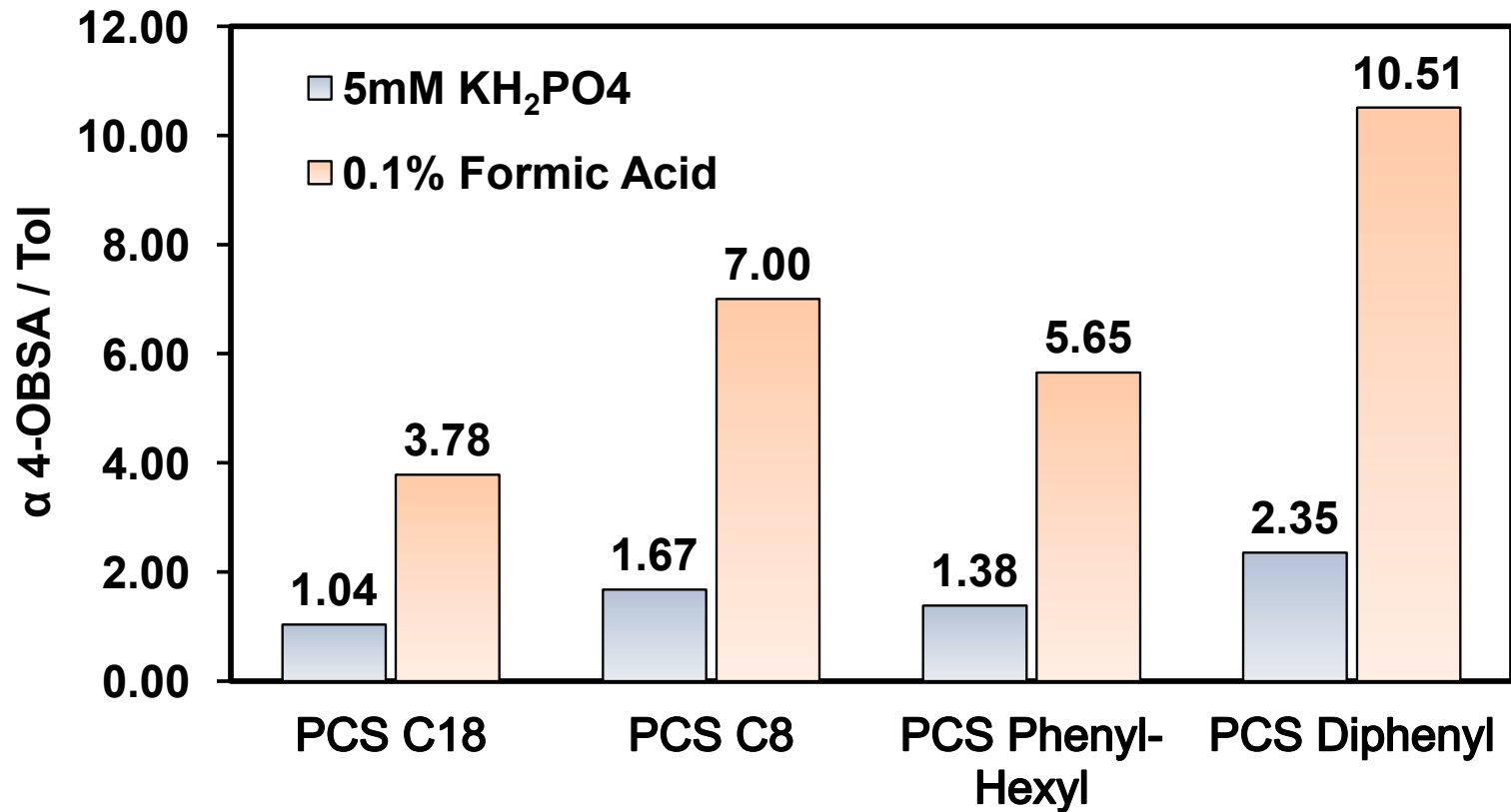

4.6x100 mm, 1.80 mL/min, 40°C, 1.0 µL Inj, 215 nm
MP – 50/50 ACN/H₂O + 5mM KH₂PO₄ pH = 2.5 or 0.1% formic acid pH = 2.7

Ionized Solute Overload Studies

4.6x100 mm, 1.20 mL/min, 30° C, 4.0 μ L Inj, 254 nm
MP - ACN/H₂O + 0.1% formic acid, Amitriptyline 3 to 10,000 ng

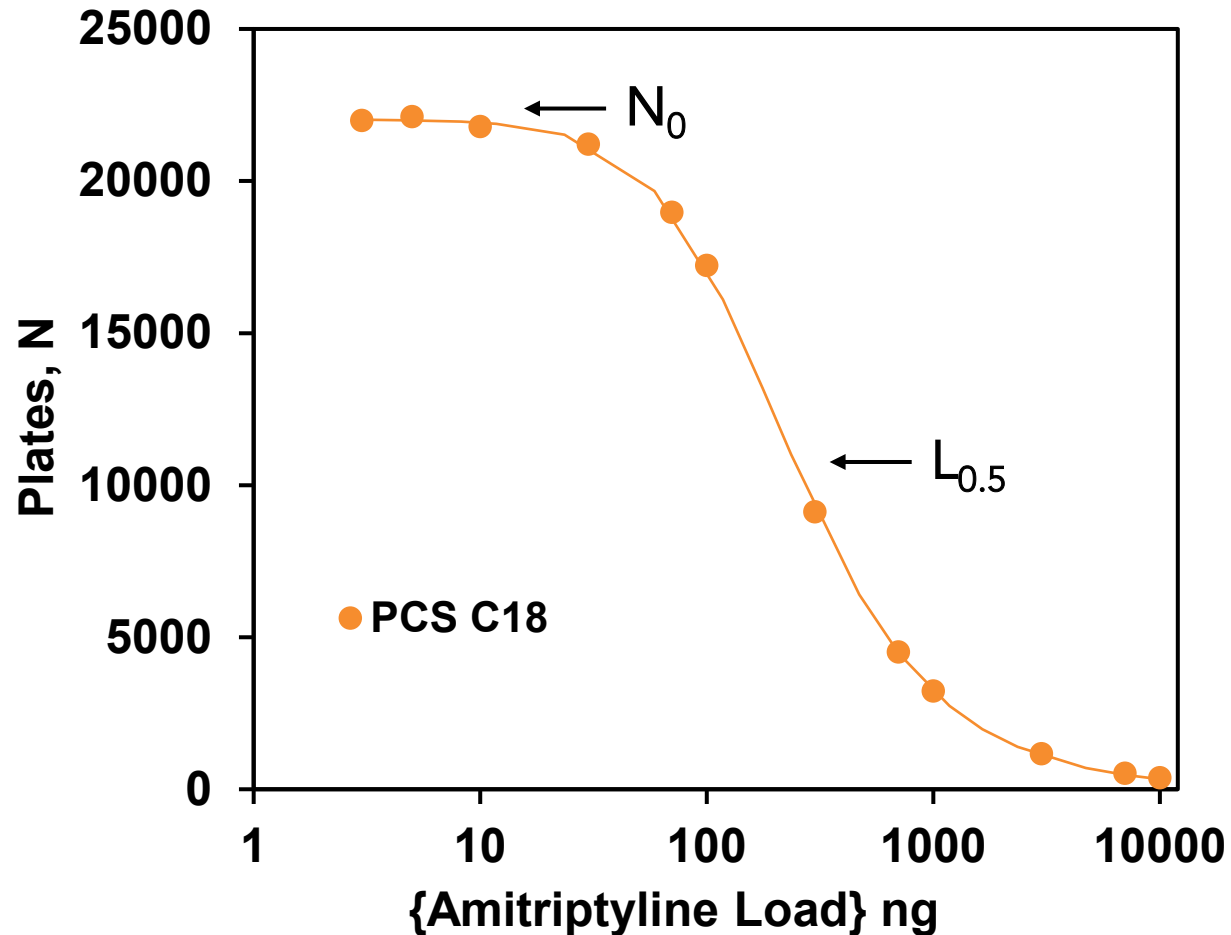

Phase	ACN %	Calculated	
		N_0	$\{L_{0.5}\}$ ng
PCS C18	19	22025	236

Evaluate ionized solute overload through [2]


- Limiting plate count (N_0)
- Load capacity at 50% of N_0 ($L_{0.5}$)

Experimental results fit to universal overload equation to solve for N_0 and $L_{0.5}$ [2]

[2] J. Dai, P.W. Carr, D.V. McCalley, A new approach to the determination of column overload characteristics in reversed-phase liquid chromatography, *Journal of Chromatography A* 1216 (2009) 2474–2482.

Ionized Solute Overload Studies

4.6x100 mm, 1.20 mL/min, 30° C, 4.0 µL Inj, 254 nm
MP - ACN/H₂O + 0.1% formic acid, Amitriptyline 3 to 10,000 ng

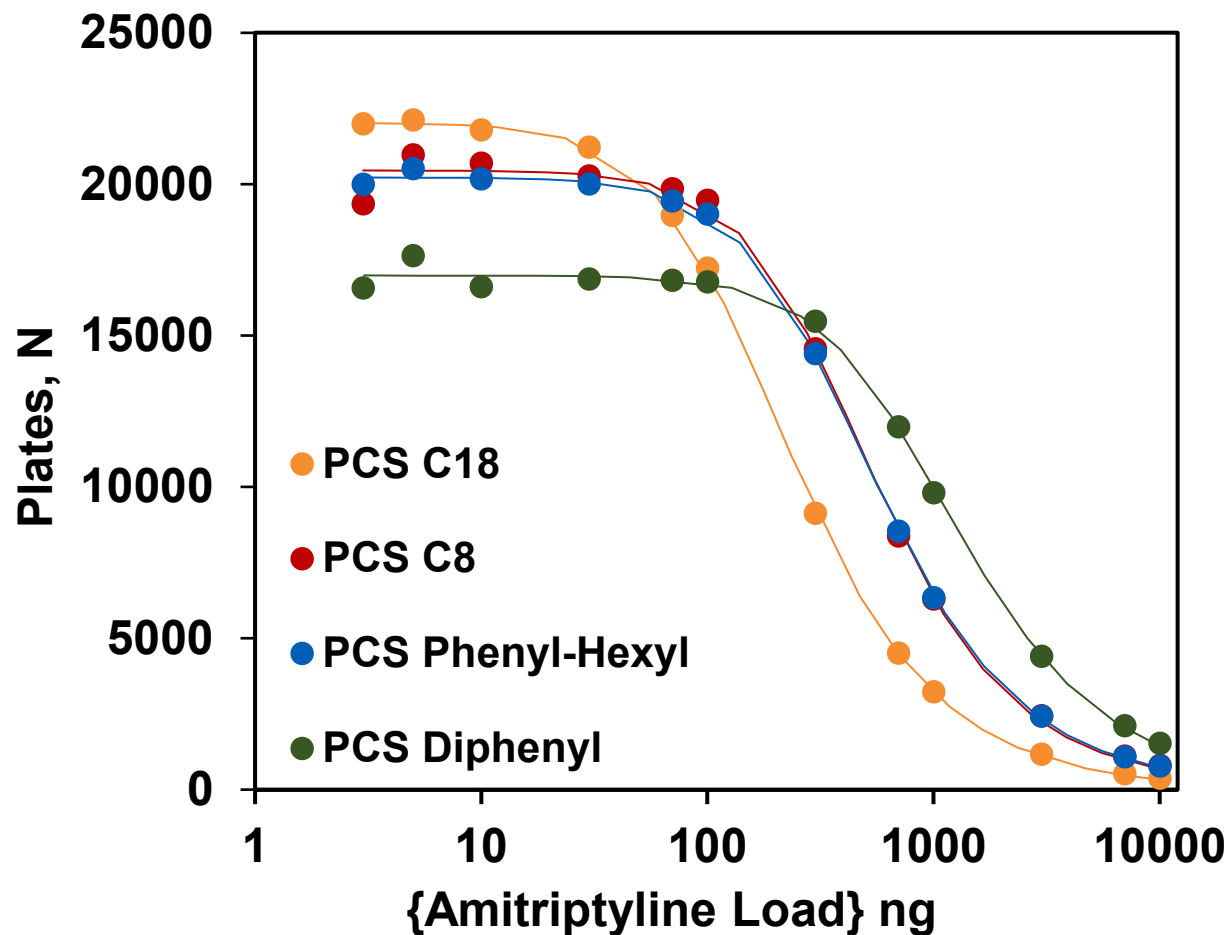

Phase	ACN	Calculated	
	%	N ₀	{L _{0.5} } ng
PCS C18	19	22025	236
PCS C8	16	20453	551
PCS Phenyl-Hexyl	16	20219	558
PCS Diphenyl	12	16984	1297

Similar trends for k' and TF

Ionized Solute Overload Studies

	PCS			
	C18	C8	Phenyl-Hexyl	Diphenyl
α 4-OBSA / Tol	3.78	7.00	5.65	10.51
Amitriptyline (+)				
N ₀	22025	20453	20219	16984
L _{0.5}	236	551	558	1297
Tetracaine (+)				
N ₀	21732	20112	18709	13476
L _{0.5}	355	868	988	1691
Naphthalene sulfonic acid (-)				
N ₀	14189	13504	14778	13961
L _{0.5}	240	483	419	769

4.6x100 mm, 1.20 mL/min, 30° C, 4.0 μ L Inj, 254 nm
MP - ACN/H₂O + 0.1% formic acid,


General trends of $\alpha_{4\text{-OBSA/Tol}}$

- Retention of ionized solutes
- Good correlation to L_{0.5} (R² = 0.9496, 0.9168, and 0.9934)

Similar basic solute results for PCS C8 and Phenyl-Hexyl, despite difference in positive character

Beta Blocker Separation

2.1x100 mm, 0.40 mL/min, 30°C, 1.0 µL Inj, 0 – 50% B in 10 min, MP A – H₂O + 0.1% FA, MP B – MeOH + 0.1%FA, LC-MS ESI conditions [3]

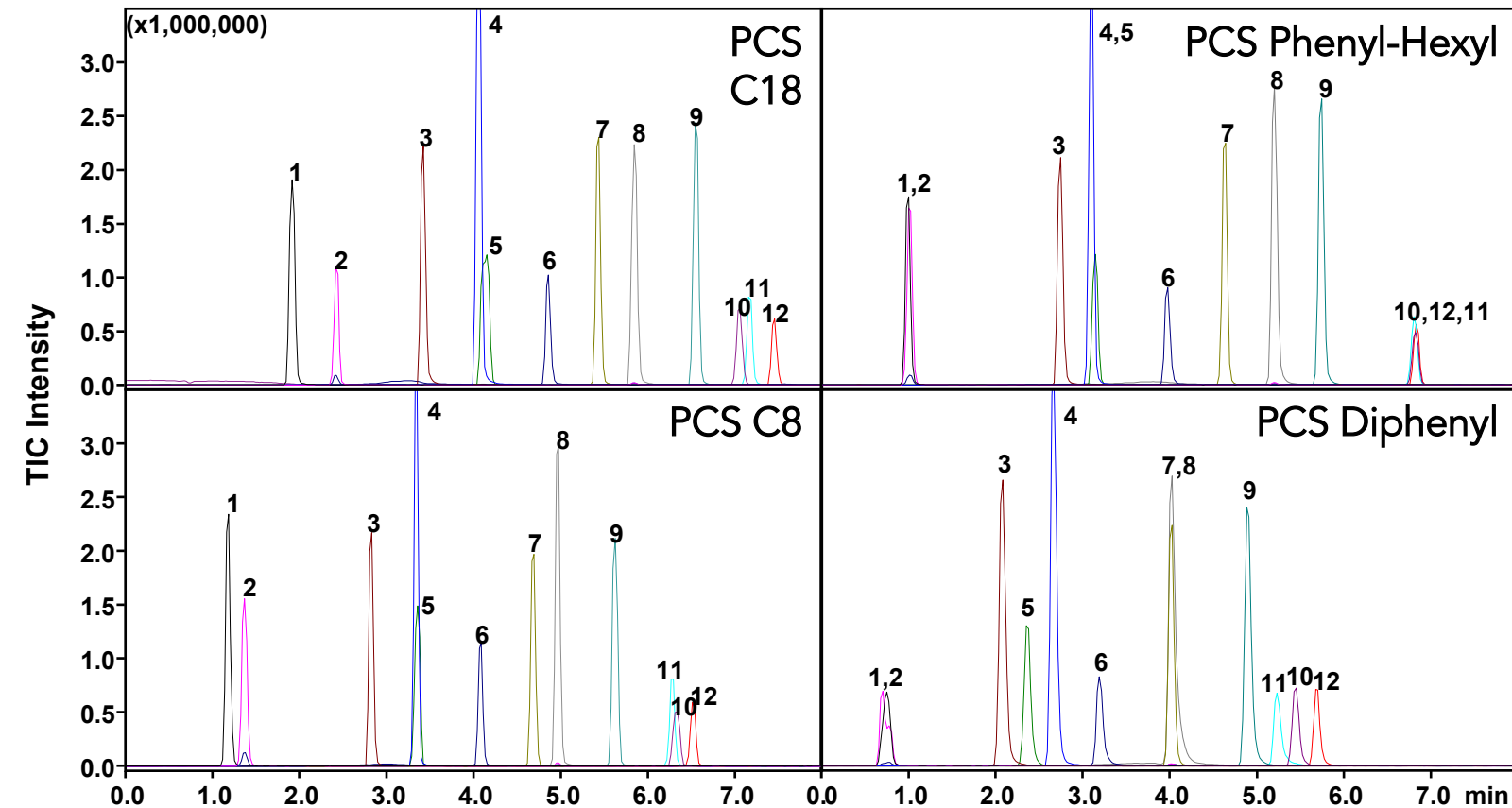

1. Sotalol 2. Atenolol 3. Pindolol 4. Carteolol 5. Nadolol 6. Metoprolol 7. Acebutolol
8. Oxprenolol 9. Bisoprolol 10. Labetalol 11. Propranolol 12. Betaxolol 0.10 ng/mL

Average retention time


- C18 > Phenyl-Hexyl > C8 > Diphenyl
- Trends with positive character
 $\alpha_{4\text{-OBSA/Tol}}$

Hydrophilic vs hydrophobic solute retention

[3] P. Szpot, K. Tusiewicz, O. Wachelko, M. Zawadzki, Application of UHPLC-QqQ-MS/MS Method for Quantification of Beta-Adrenergic Blocking Agents (β -Blockers) in Human Postmortem Specimens, *Molecules* 29 (2024) 4585.

Beta Blocker Separation

2.1x100 mm, 0.40 mL/min, 30°C, 1.0 µL Inj, 0 – 50% B in 10 min, MP A – H₂O + 0.1% FA, MP B – MeOH + 0.1%FA, LC-MS ESI conditions [3]

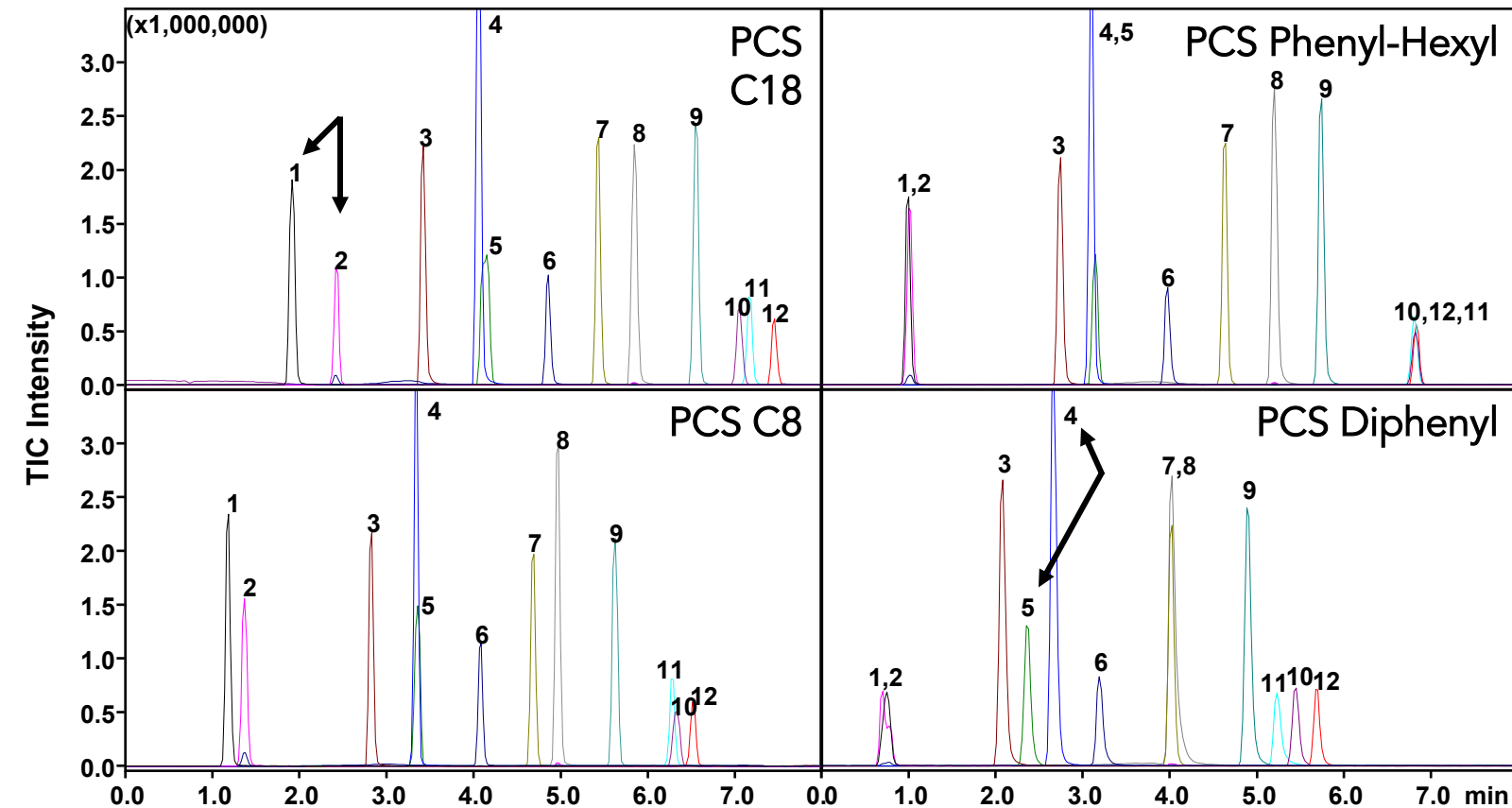

1. Sotalol 2. Atenolol 3. Pindolol 4. Carteolol 5. Nadolol 6. Metoprolol 7. Acebutolol
8. Oxprenolol 9. Bisoprolol 10. Labetalol 11. Propranolol 12. Betaxolol 0.10 ng/mL

Changes in selectivity observed

- Importance of primary hydrophobic ligand

Peaks 1 and 2 PCS C18

Peaks 4 and 5 PCS Diphenyl

[3] P. Szpot, K. Tusiewicz, O. Wachelko, M. Zawadzki, *Application of UHPLC-QqQ-MS/MS Method for Quantification of Beta-Adrenergic Blocking Agents (β -Blockers) in Human Postmortem Specimens, Molecules* 29 (2024) 4585.

Conclusions

Primary hydrophobic ligand does affect the positive character of these cation modified stationary phases.

- Controlling for surface area and concentration of the positive charge

$\alpha_{4\text{-OBSA/Tol}}$ classifies HALO PCS columns as low to medium positive character, as designed [1].

As positive character of the phase increases, an increase in resistance to the effects of ionized solute overloading ($L_{0.5}$, tailing factor, and k') was observed.

Anionic exchange parameter was found to correlate to general retention and selectivity trends of ionized acids and bases.

Acknowledgments

Research and Development Team


- Mark Haynes
- Bryant McKnight
- Timothy Langlois

Engineering Team

- Matt Jackson

Thank you for your time today!

Questions

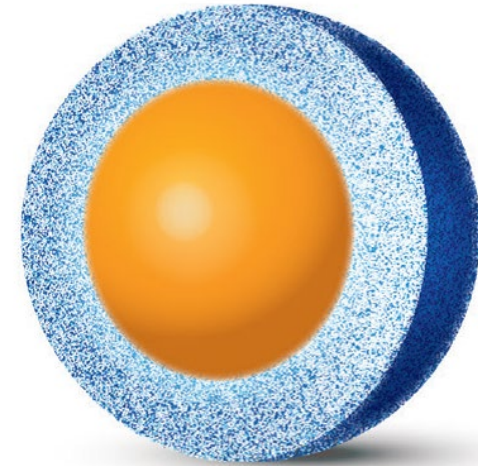



advancedmaterialstechnology



halocolumns.com



3521 Silverside Road, Suite 1-K
Quillen Building
Wilmington, DE 19810




(302) 992-8060



Ionized Solute Overload Studies

4.6x100 mm, 1.20 mL/min, 30° C, 4.0 µL Inj, 254 nm
MP - ACN/H₂O + 0.1% formic acid, Amitriptyline 3 to 10,000 ng

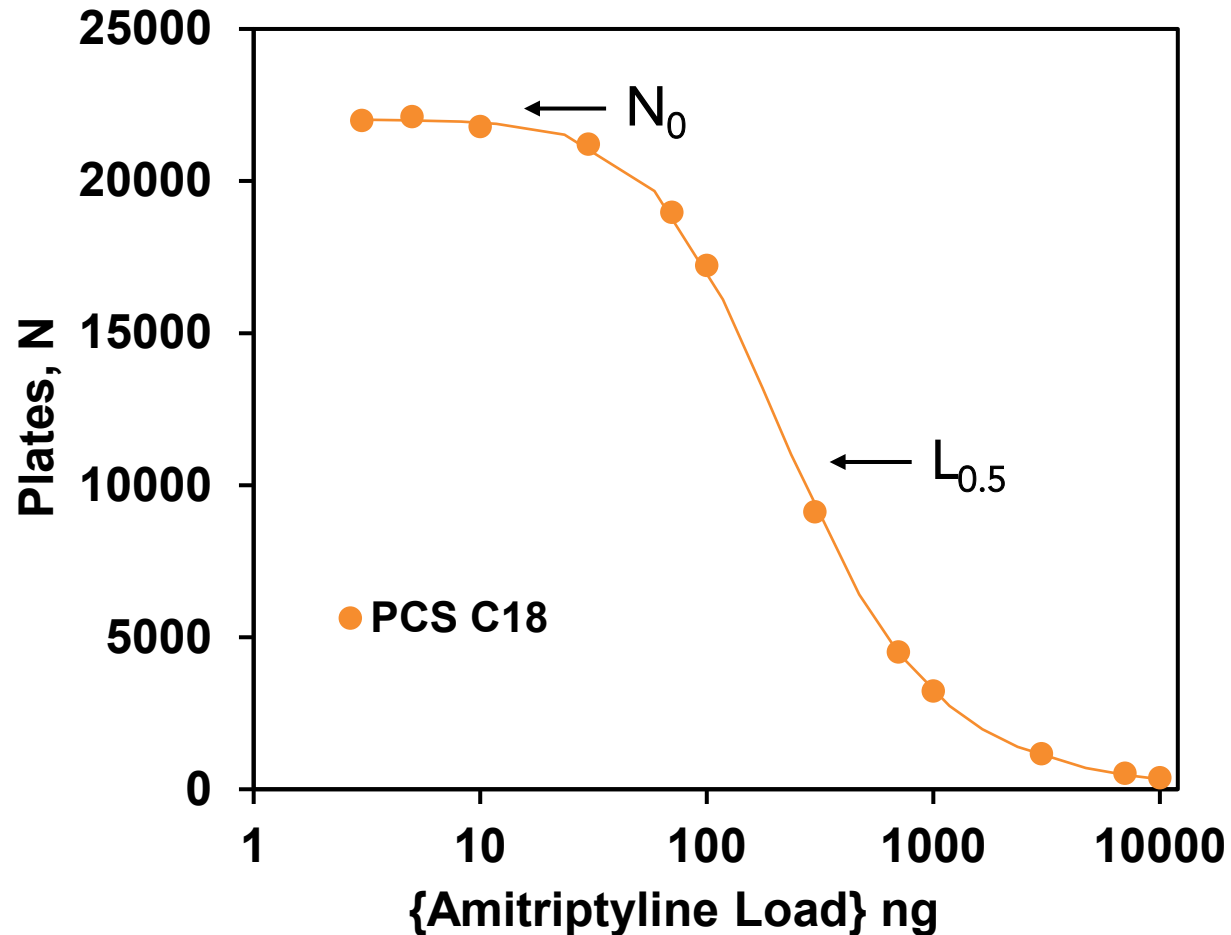

Phase	ACN %	Calculated	
		N_0	$\{L_{0.5}\}$ ng
PCS C18	19	22025	236

Experimental results fit to universal overload equation to solve for N_0 and $L_{0.5}$ [2]

$$\frac{N}{N_0} = \frac{1 + bL'}{1 + bL' + (b + 1)L'^2}$$


[2] J. Dai, P.W. Carr, D.V. McCalley, A new approach to the determination of column overload characteristics in reversed-phase liquid chromatography, *Journal of Chromatography A* 1216 (2009) 2474–2482.

Anionic Exchange Measurement

4.6x100 mm, 1.80 mL/min, 40° C, 1.0 µL Inj, 215 nm
MP – 50/50 ACN/H2O + 5mM KPO4 pH = 2.5

1.Toluene (0.50 mg/mL)

2. 4-OBSA (0.50 mg/mL)

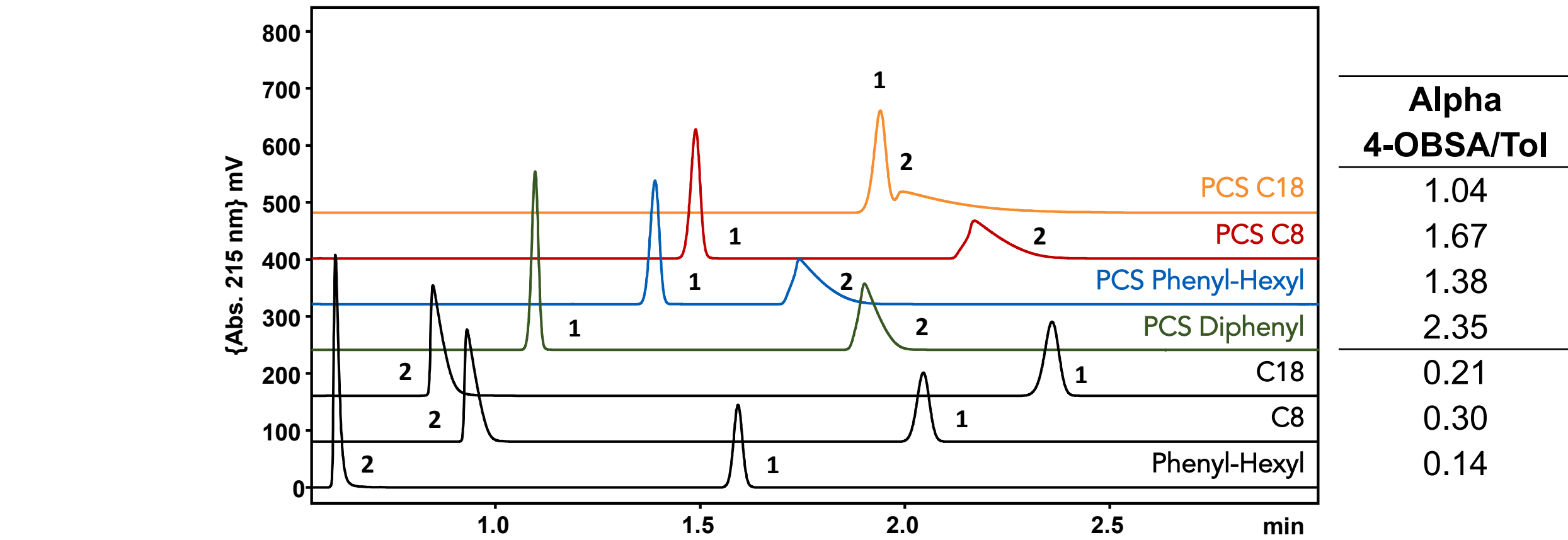


Anionic Exchange Measurement

4.6x100 mm, 1.80 mL/min, 40° C, 1.0 μ L Inj, 215 nm
MP – 50/50 ACN/H₂O + 5mM KH₂PO₄ pH = 2.5 , Ref [1]


1.Toluene (0.50 mg/mL)

2. 4-OBSA (0.50 mg/mL)

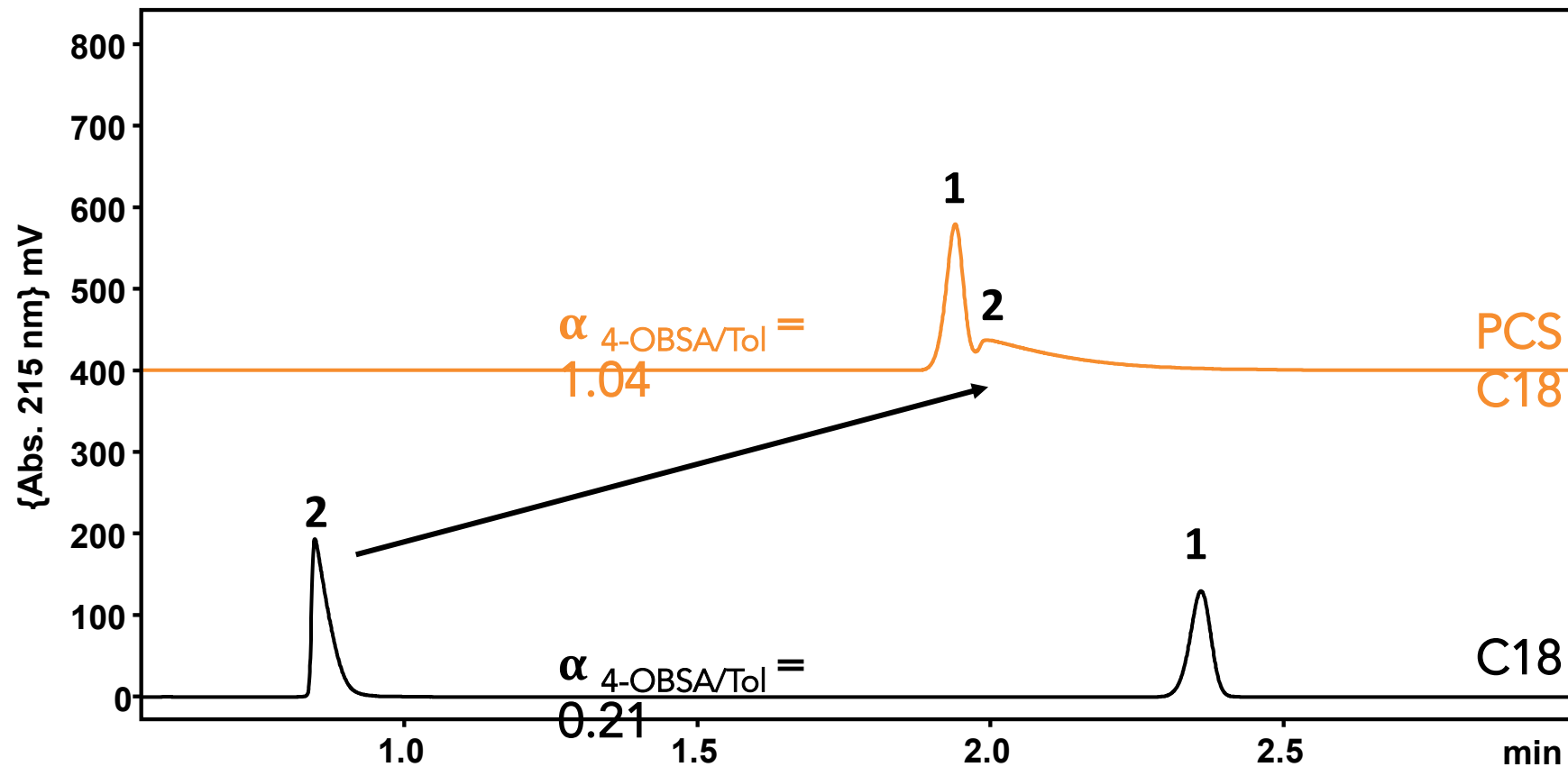

Ionized Basic Solute Overload

4.6x100 mm, 1.20 mL/min, 30° C, 4.0 µL Inj, 254 nm
MP - ACN/H₂O + 0.1 % formic acid (v/v),
Amitriptyline 3 to 100 ng

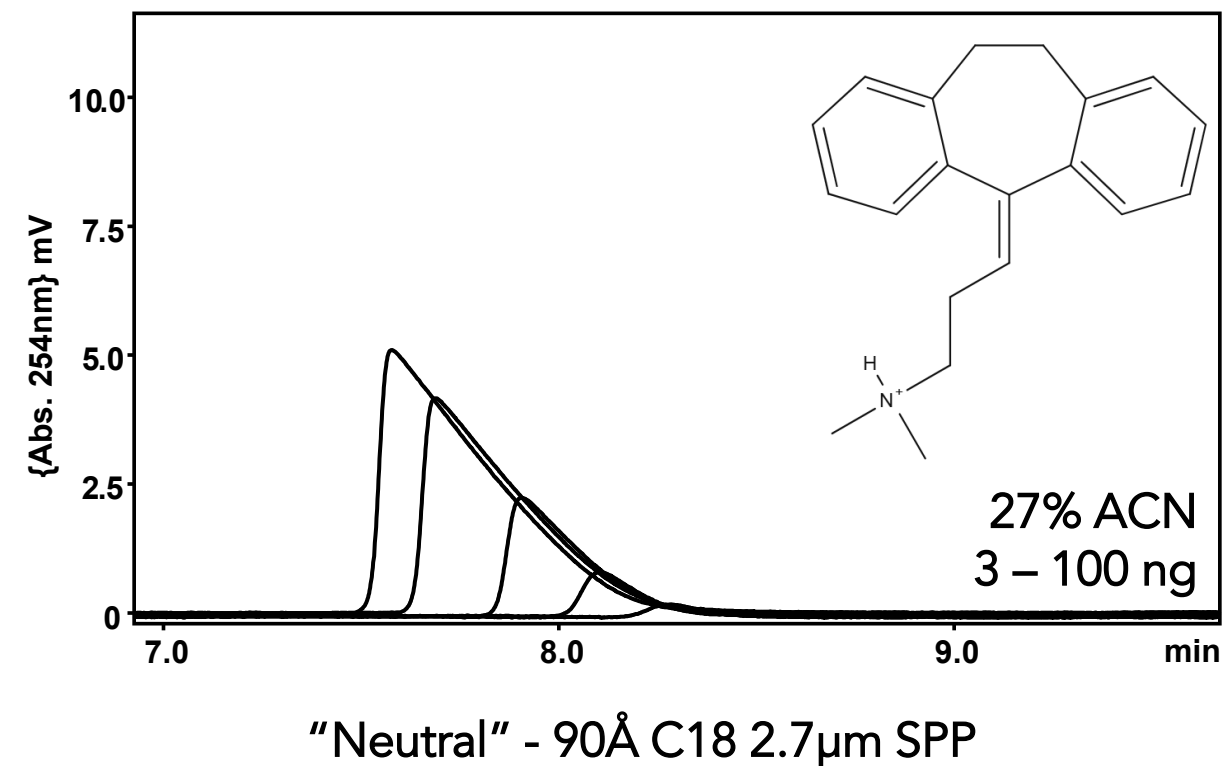

Overload [Base] <<<< [Neutral]

Non-linear chromatography

Tailing Factor ↑ (1.50 → 5.43)

Plates, N ↓ (17,610 → 3,423)

Retention Factor, k' ↓ (11.1 → 10.1)